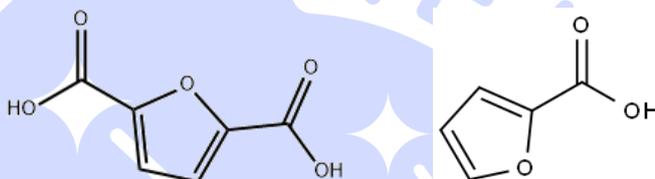
Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Furandicarboxylic acid 呋喃二甲酸	Sample properties 样品性状	Liquid 固体
Date of receipt 收样日期	2025/09/02	During testing 测试期间	2025/09/07 ~ 2025/09/08
Test the ingredients and structural formula			
测试成分及结构式			
 <p>Furandicarboxylic acid 呋喃二甲酸</p> <p>Furoic acid 呋喃甲酸</p>			
Experimental Requirements			
实验要求			
To develop a method for determining furandicarboxylic acid content with separation from its main impurity (furoic acid) and a resolution greater than 1.5 开发方法测定呋喃二甲酸含量, 要求与主要杂质呋喃甲酸分离, 分离度大于 1.5			
Reference Method			
参考方法			
/			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
Water 水	Ultrapure water 超纯水	Welch 月旭	

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile phase A: Weigh 2.72 g of potassium dihydrogen phosphate, dissolve it in 1000 ml of pure water, adjust the pH to 2.2 with phosphoric acid after full dissolution, mix well and filter to obtain the solution.

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Pure water.

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1.2.3. Test solution (供试品溶液):

Test solution L or S: Weigh 10 mg of the test sample L or S into a 10 ml volumetric flask, dissolve it with pure water, dilute to the marked line, and mix well to obtain.

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1.2.4. Furoic acid reference solution: Take 10 mg of the reference substance into a 10 ml volumetric flask, dissolve it with pure water, dilute to the marked line, and mix well to obtain.

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Tel: 400-810-6969

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Web: www.welchmat.com

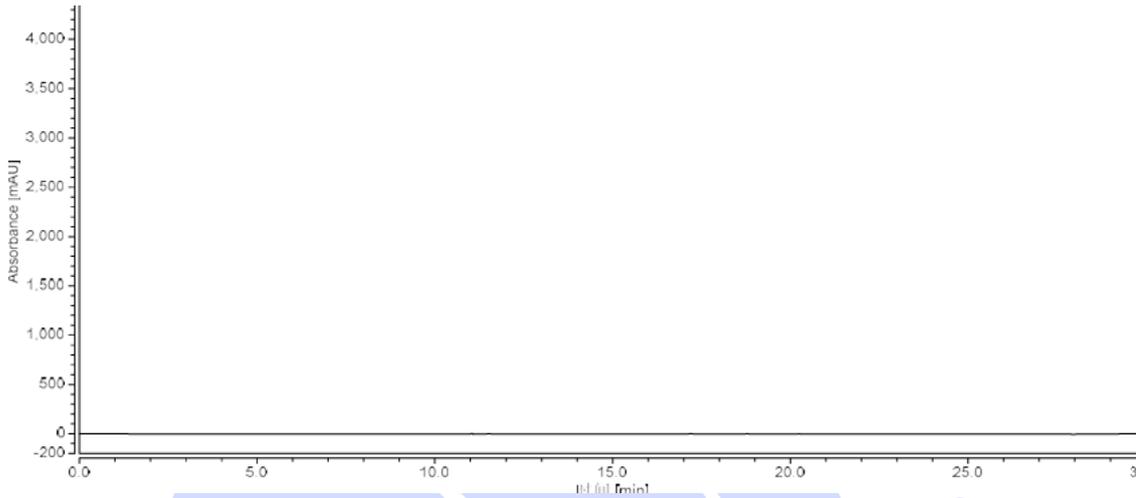
1.2.5. Mixed solution: Take equal proportions of the test solution and the reference solution of the impurity furancarboxylic acid, mix them well to obtain.

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
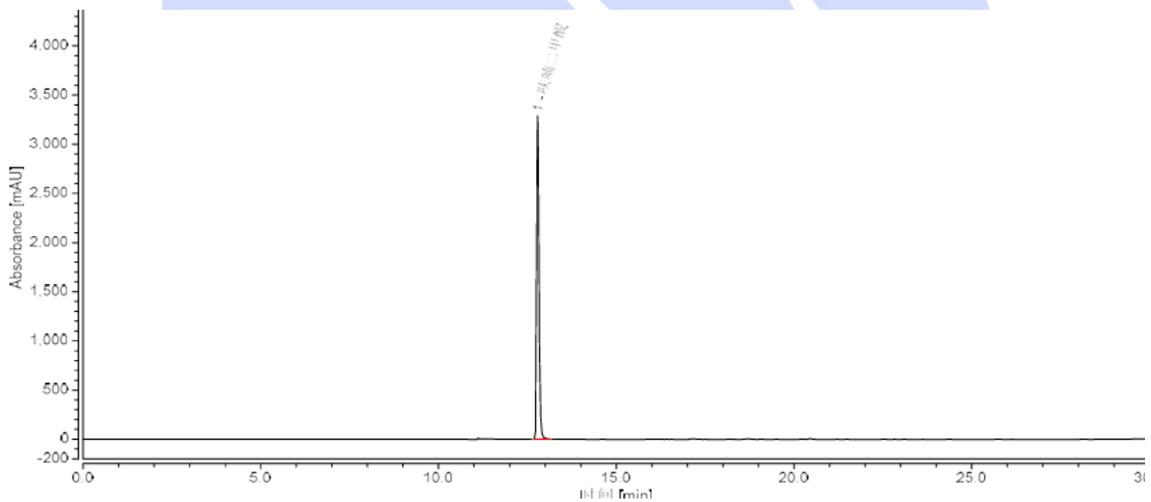
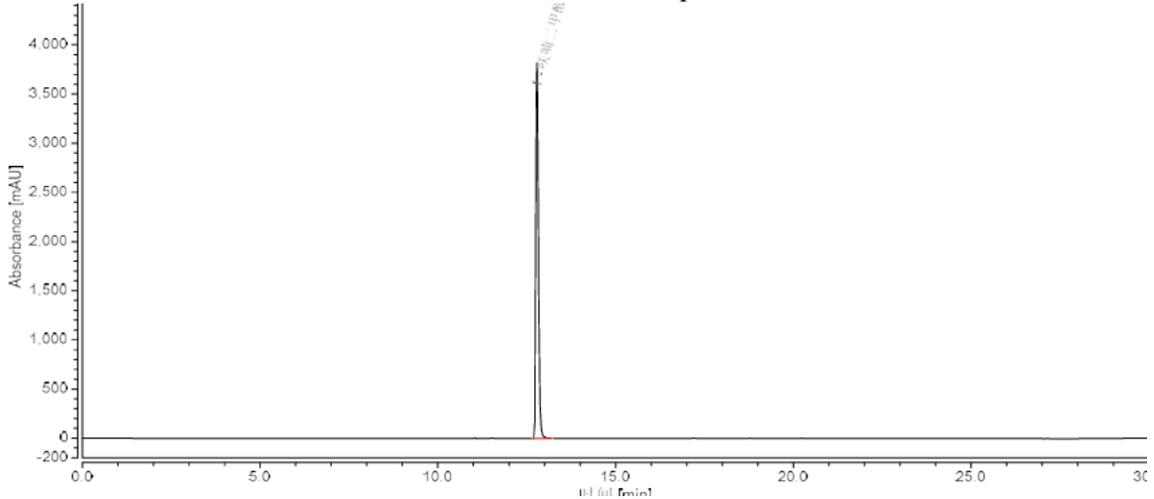
(3) Analysis spectrum of Furoic acid reference solution:

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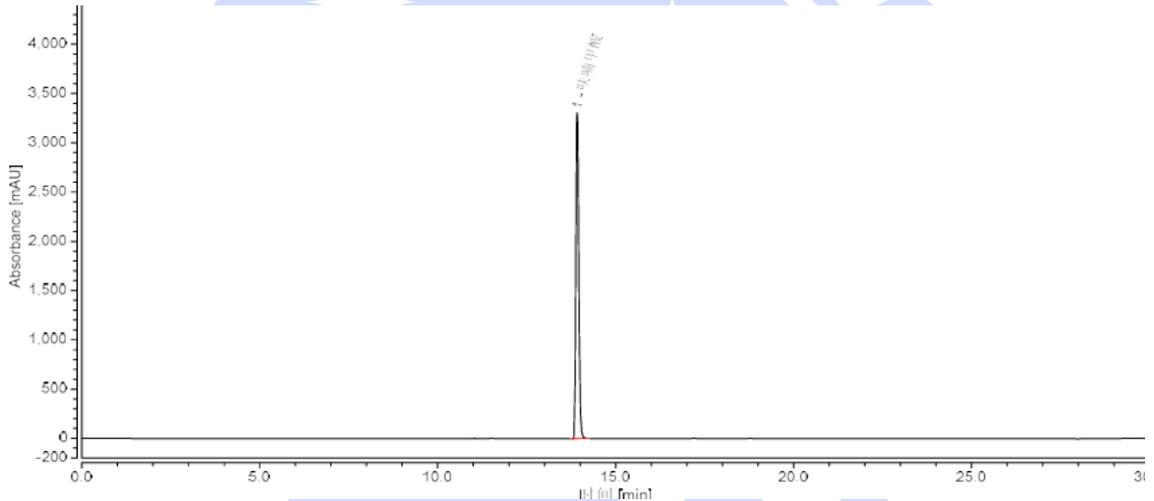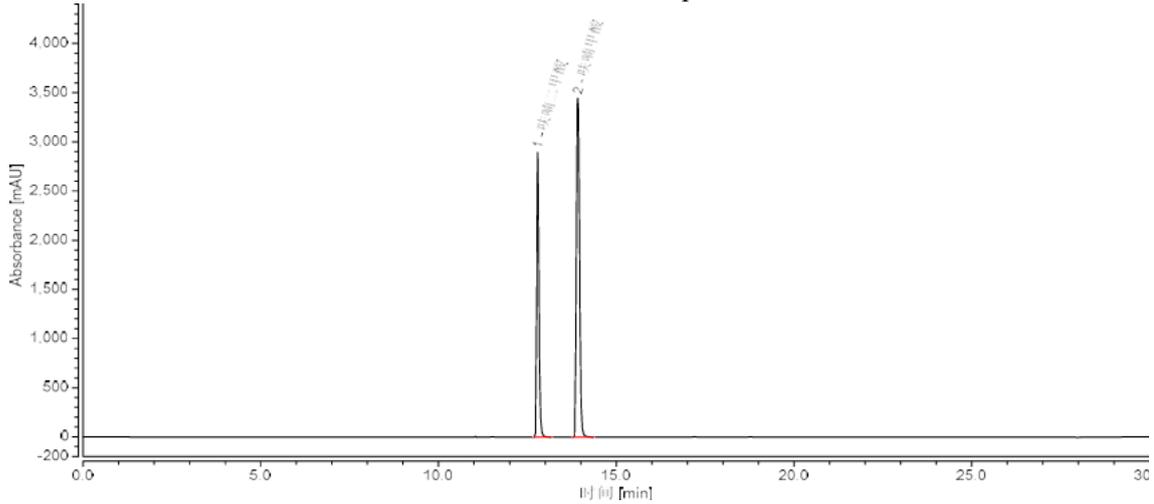
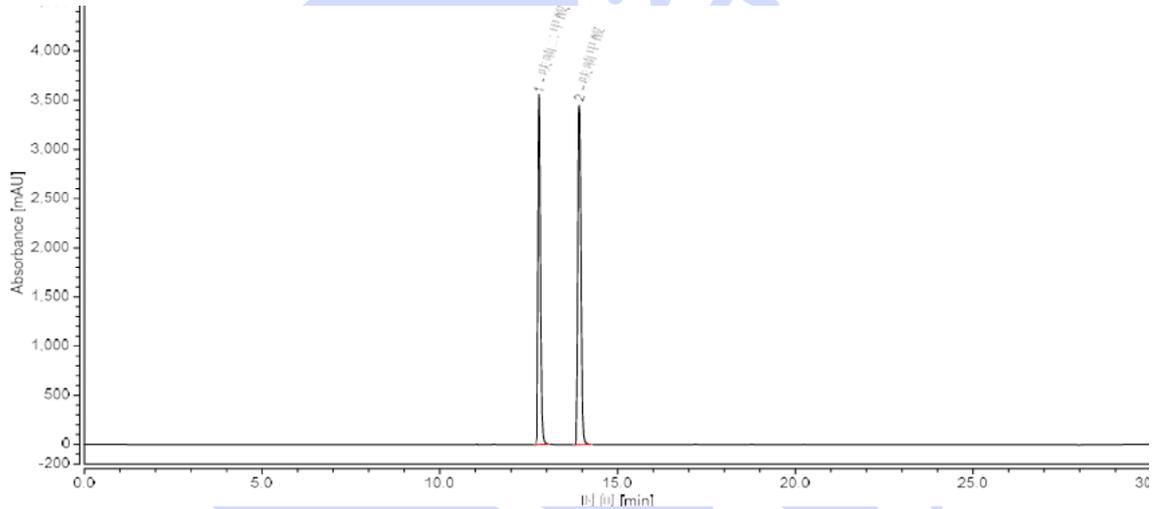
3. Conclusion (结论)

Using Welch Ultimate LP-C18 (4.6×250mm, 5µm) analyzed the content of Furandicarboxylic acid under this chromatographic condition. The resolution between the target compound and the impurity Furoic acid is better than 7.0 , which meets the customer's testing requirements.

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