Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Protein Content 蛋白含量	Sample properties 样品性状	Solid 固体
Date of receipt 收样日期	2025/09/25	During testing 测试期间	2025/09/29 ~ 2025/10/09
Test the ingredients and structural formula			
测试成分及结构式			
/			
Experimental Requirements			
实验要求			
Develop a method to analyze protein content in samples using bovine serum albumin (BSA) as a reference. 开发方法, 以牛血清白蛋白为参照测定样品中蛋白含量			
Reference Method			
参考方法			
/			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
Water 水	Ultrapure water 超纯水	Welch 月旭	
Acetonitrile 乙腈	HPLC 色谱级	Welch 月旭	
Potassium Dihydrogen Phosphate	AR 分析级	Aladdin 阿拉丁	

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile Phase: Weigh 13.6 g of potassium dihydrogen phosphate into 1000 mL of pure water, stir

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Tel:400-810-6969

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Web:www.welchmat.com

until fully dissolved. Adjust the pH to 6.8 using a potassium hydroxide solution, mix well and then obtain the mobile phase by suction filtration.

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1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

Acetonitrile (乙腈); Mobile Phase used for dilution (稀释时可以用流动相)

1.2.3. BSA reference solution: Weigh 10 mg of the reference standard into a 100 ml volumetric flask, dissolve it with mobile phase, dilute to the marked volume, and mix well to obtain the solution.

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1.2.4. BSA Sensitivity Solution: Pipette 0.5 mL of the reference solution into a 100 mL volumetric flask, dilute to the marked volume with the mobile phase, and mix well to obtain the solution.

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1.2.5. Sample Solutions: Weigh 10 mg each of Sample 1, Sample 2, and Enzyme Sample separately into 10 mL volumetric flasks, dilute to the marked volume with a solvent, and mix well to obtain the solutions.

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1.2.6. Spiked Solutions: Weigh Sample 1 and Sample 2 separately, prepare them into solutions with a concentration of 20 mg/mL, and then add an appropriate amount of protein reference solution to these solutions to obtain spiked solutions with a protein concentration of 5 µg/mL.

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2. Chromatogram and data (谱图和数据)

(1) Analysis spectrum of Blank solvent solution (空白溶剂分析图谱)

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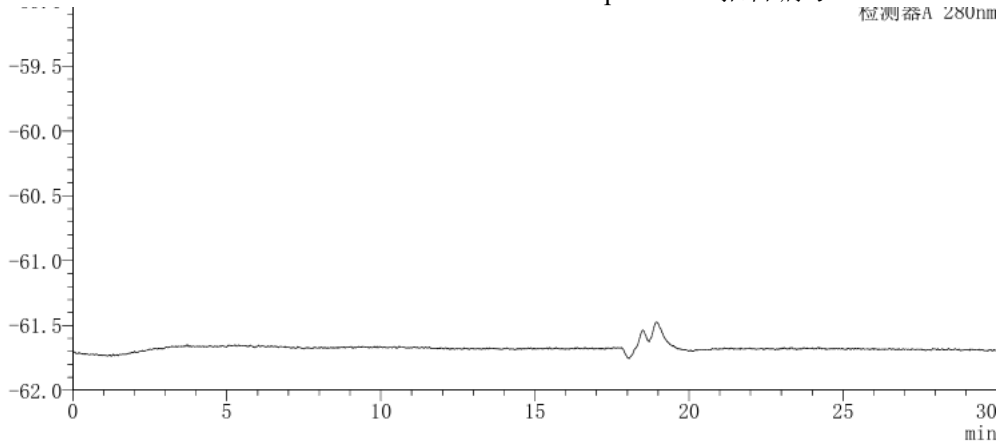
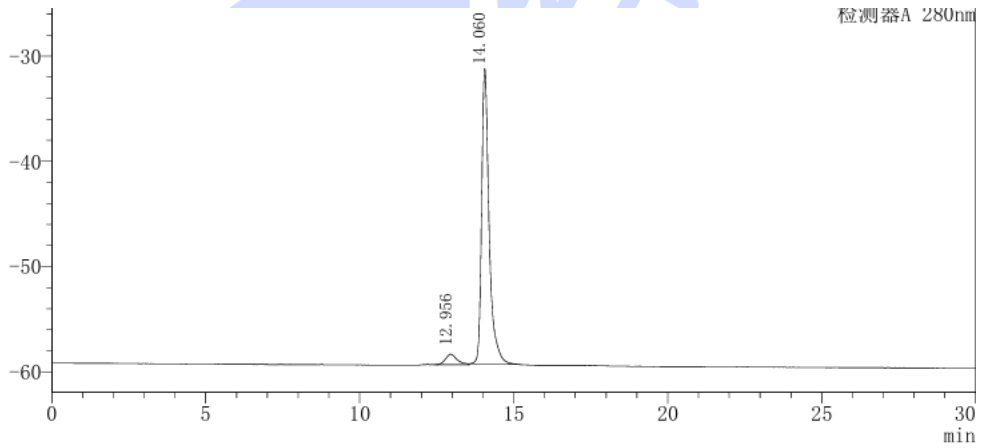
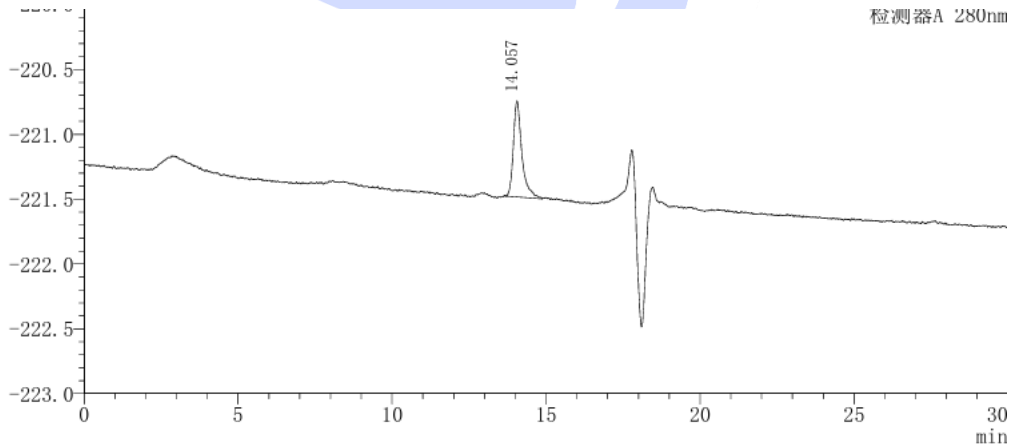
(2) Analysis spectrum of BSA Sensitivity solution (牛血清白蛋白灵敏度溶液分析图谱):

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
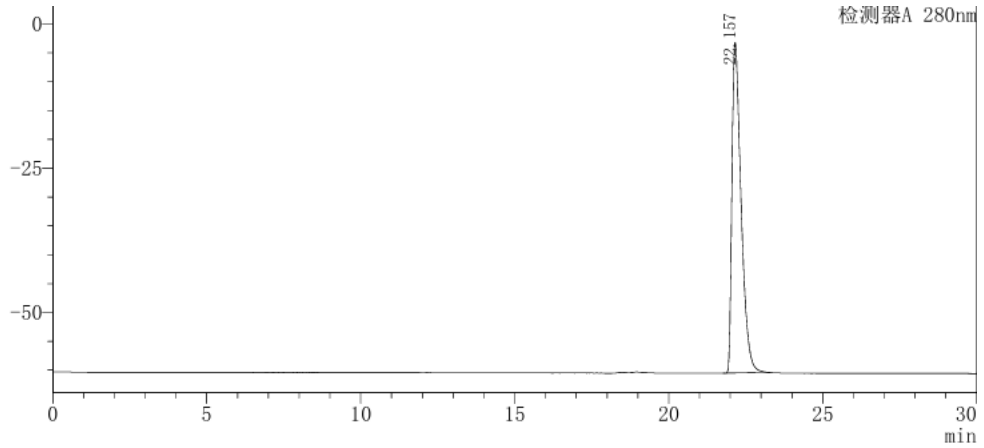
(3) Analysis spectrum of Sample 1 solution (样品 1 溶液分析图谱):

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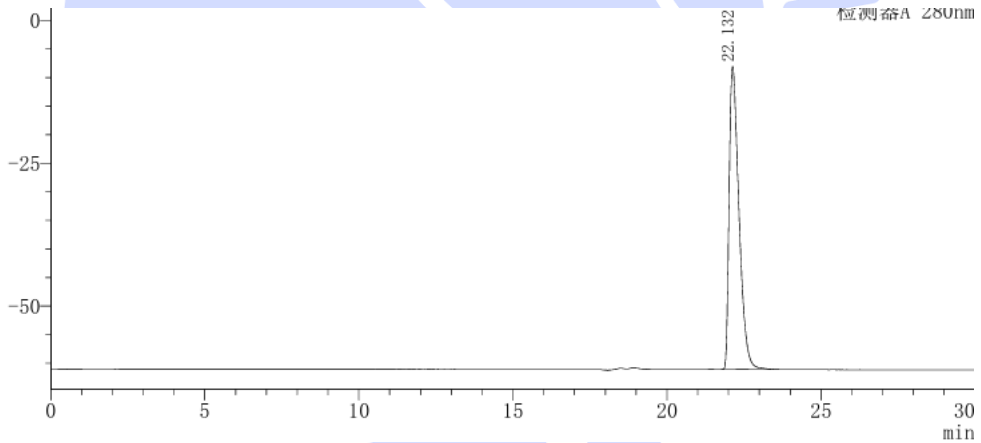(4) Analysis spectrum of Sample 2 solution (样品 2 溶液分析图谱):

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(5) Analysis spectrum of Enzyme Sample solution (酶样品溶液分析图谱):

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
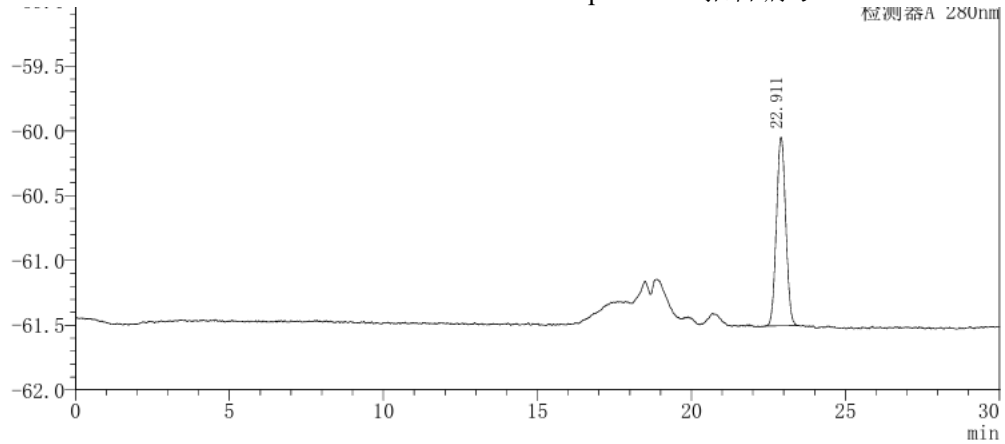
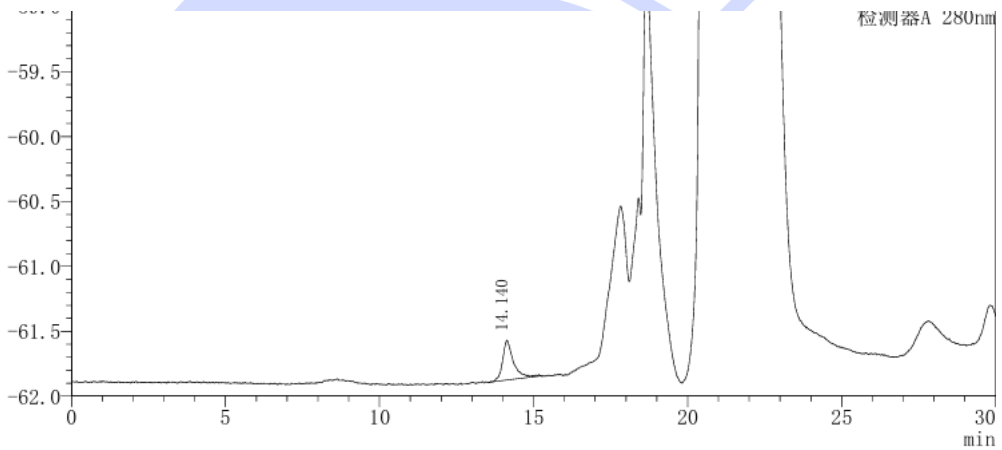
(6) Analysis spectrum of Sample 1 spike solution (样品1加标溶液分析图谱):

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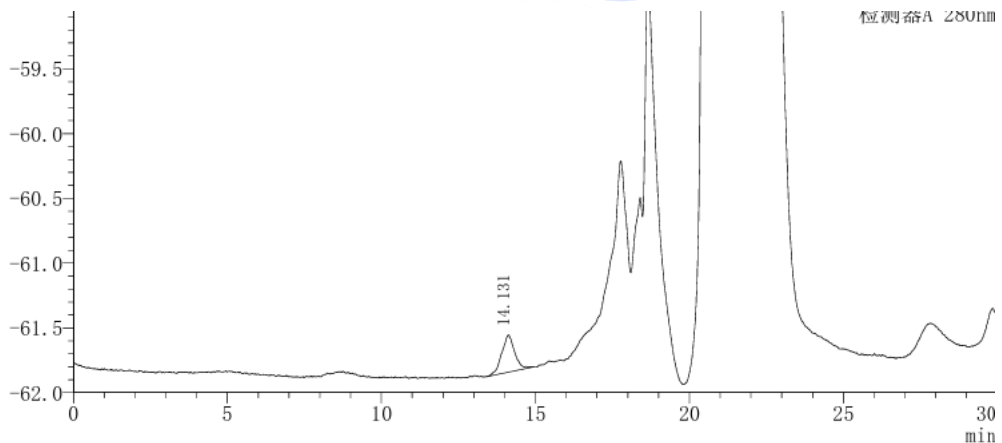(7) Analysis spectrum of Sample 2 spike solution (样品2加标溶液分析图谱):

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3. Conclusion (结论)

Using Welch Xtimate Bio SEC-300 (7.8×0mm, 3μm) analyzed the content of Protein under this chromatographic condition , the result meets the customer's testing requirements.

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