Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Aminophenazone, benzoic acid, sorbic acid, saccharin and dehydroacetic acid 安赛蜜、苯甲酸、山梨酸、糖精钠 和脱氢乙酸	Sample properties 样品性状	Colorless transparent liquid无色透明液体
Date of receipt 收样日期	2025/09/04	During testing 测试期间	2025/09/04 ~ 09/09
Test the ingredients and structural formula			
测试成分及结构式			
Aminophenazone, benzoic acid, sorbic acid, saccharin and dehydroacetic acid 安赛蜜、苯甲酸、山梨酸、糖精钠和脱氢乙酸			
Experimental Requirements			
实验要求			
Separation meets requirements. 分离满足要求。			
Reference Method			
参考方法			
GB5009.28-2016			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
乙酸铵	AR	阿拉丁	
乙酸	AR	国药	

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

20mM ammonium acetate solution: 1.54g of ammonium acetate was weighed, 1000ml of water was added to dissolve it, and the pH value was adjusted to 6.7 with acetic acid.

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1.2.2. Test solution (供试品溶液):

We accurately weighed approximately 2 g of the sample and transferred it to a 50 mL screw-capped centrifuge tube. After adding about 25 mL of water and vortex-mixing, the sample was subjected to 50°C water bath ultrasonication for 20 minutes. Following cooling to room temperature, 2 mL of potassium ferrocyanide solution and 2 mL of zinc acetate solution were added. The mixture was thoroughly mixed and centrifuged at 8000 r/min for 5 minutes. The aqueous phase was then transferred to a 50 mL volumetric flask. Adding 20 mL of water to the residue, vortex-mixing followed by 5 minutes of ultrasonication was performed, after which the flask was centrifuged at 8000 r/min for 5 minutes. The aqueous phase was again transferred to the same 50 mL volumetric flask and made up to the mark with water. A portion of the supernatant was filtered through a 0.22 μ m membrane for liquid chromatography analysis..

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1.2.3. Reference solution (对照溶液):

Mix of asemit, benzoic acid, sorbic acid, and saccharin sodium: provided by the customer at a concentration of 100 μ g/mL.

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Dehydroacetic acid: supplied by the customer, concentration of 100 μ g/mL

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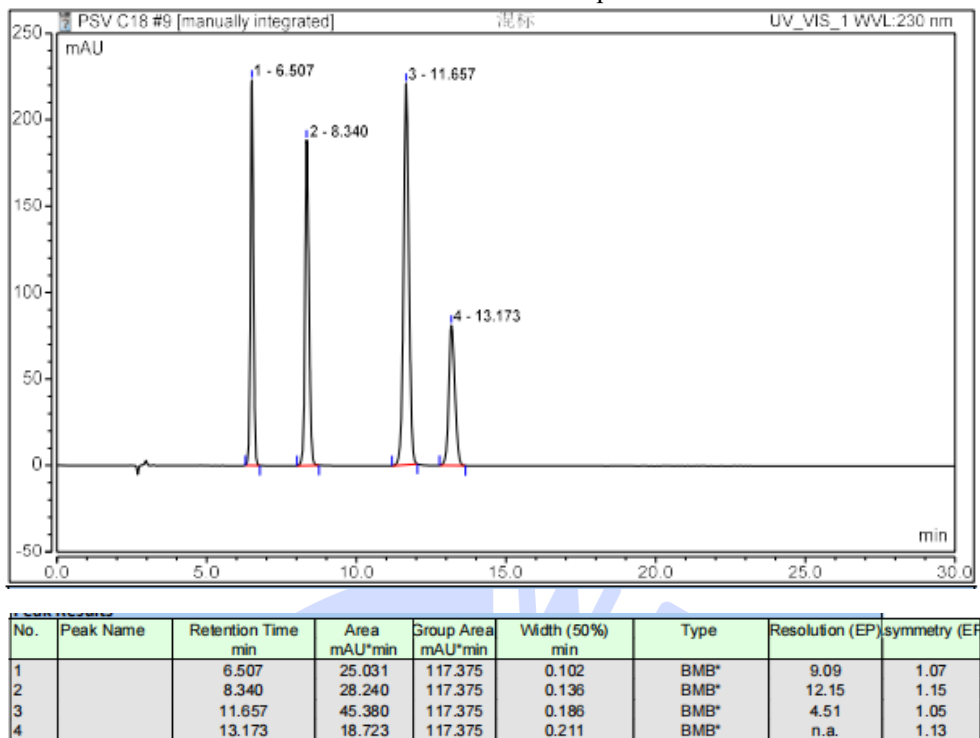
2. Chromatogram and data (谱图和数据)

(1) Reference solution (对照溶液图)

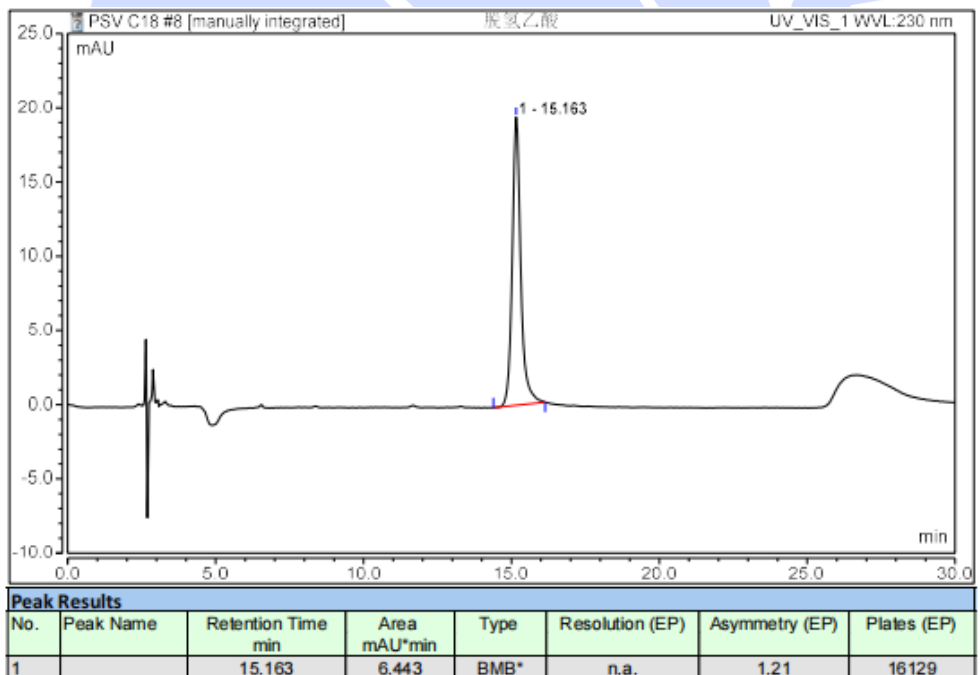
Spectra of mixtures of amines, benzoic acid, sorbic acid and sodium saccharin.

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Dehydroacetic acid solution diagram

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