Test report

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Sample Information 样品信息			
Sample Name 样品名称	ANEMARRHENAE RHIZOMA 知母	Sample properties 样品性状	Presenting a light yellow elongated shape 呈淡黄色长条状
Date of receipt 收样日期	2025/09/01	During testing 测试期间	2025/09/18 ~ 09/19
Test the ingredients and structural formula 测试成分及结构式			
New mangoside, mangoside 新芒果苷、芒果苷			
Experimental Requirements 实验要求			
Theoretical plate count should be at least 6000 based on the peak of mangiferin. 理论板数按芒果苷峰计算应不低于 6000			
Reference Method 参考方法			
Chinese Pharmacopoeia 中国药典			
Reagent information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
Phosphoric acid 磷酸	85%	Aladdin 阿拉丁	
Acetonitrile	Chromatographic pure	Welch	

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Take 2mL of phosphoric acid and add it to 998mL of water. mix well, ultrasonic degassing to obtain;
取 2mL 磷酸, 加入 998mL 水, 混匀, 超声脱气即得;

1.2.2. Blank solution (空白溶液):

50% methanol.

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1.2.3. Reference solution (对照品溶液):

Take a suitable amount of new mangiferin reference substance and mangiferin reference substance, weigh accurately, and add 50% methanol to prepare a mixed solution containing 75 μ g of new mangiferin and 50 μ g of mangiferin per 1ml.

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1.2.4. Test sample solution (供试品溶液):

Take about 0.2g of the powdered product (passed through a No. 3 sieve), weigh it precisely, and place it in a stoppered conical flask. Accurately add 50 ml of 50% methanol, seal tightly, and weigh it again. Ultrasonicate for 40 minutes (power 400W, frequency 40kHz), allow to cool, and weigh again. Use 50% methanol to make up for the lost weight, shake well, filter, and collect the filtrate to obtain the solution.

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