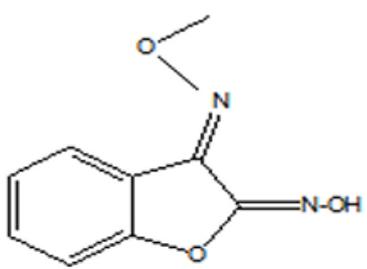
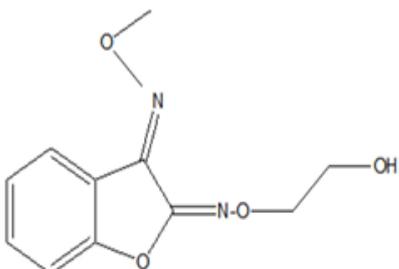
Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Benzofurans 苯并呋喃类物质	Sample properties 样品性状	Liquid 固体
Date of receipt 收样日期	2025/09/26	During testing 测试期间	2025/09/29 ~ 2025/10/09
Test the ingredients and structural formula			
测试成分及结构式			
 <p>192.17 I-4</p>		 <p>236.23 I-5</p>	
Experimental Requirements			
实验要求			
<p>Develop a method to analyze substances I-4, I-5 and their respective cis-trans isomers, with the requirements that the resolution is greater than 2 and the peak time is within 10 minutes.</p> <p>开发方法分析物质 I-4、I-5 及其各自对应的顺反异构体，要求分离度大于 2，出峰时间 10min 内</p>			
Reference Method			
参考方法			
/			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
Water	Ultrapure water	Welch	

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## 1.2. Solution preparation (溶液配制)

### 1.2.1. Preparation of mobile phase (流动相配制):

Mobile phase A: Take 1ml Phosphoric acid in 1000 ml of pure water, mix well and filter to obtain.

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### 1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

Initial proportion mobile phase (初始比例流动相)

1.2.3. I-4 reference solution: Weigh 10 mg of the reference standard into a 100 ml volumetric flask, dissolve it with a solvent, dilute to the marked volume, and mix well to obtain the solution.

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1.2.4. I-5 reference solution: Weigh 10 mg of the reference standard into a 100 ml volumetric flask, dissolve it with a solvent, dilute to the marked volume, and mix well to obtain the solution.

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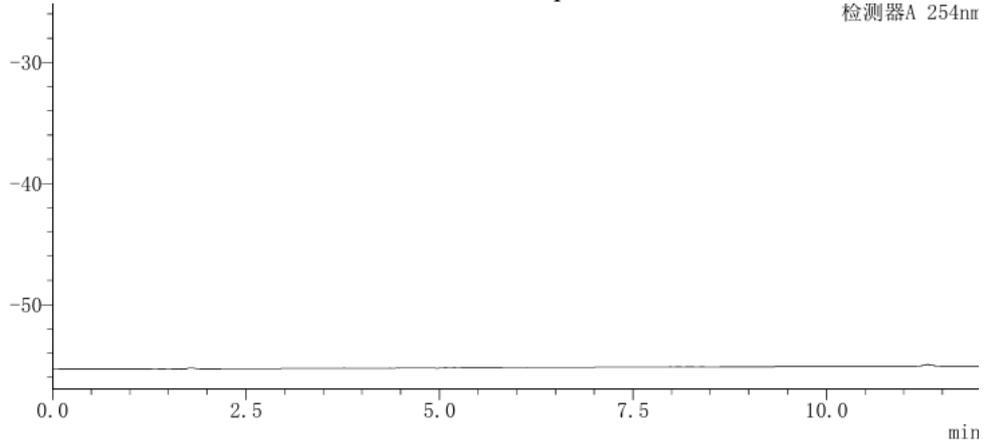
1.2.5. Mixed reference solution: Take 1 ml each of the reference solutions of I-4 and I-5 into a 10 ml volumetric flask, dilute to the marked volume with a solvent, and mix well to obtain the solution..

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
## 2. Chromatogram and data (谱图和数据)

### (1) Analysis spectrum of Blank solvent solution (空白溶剂分析图谱)

## (2) Analysis spectrum of I-4 reference solution (I-4 对照溶液分析图谱):

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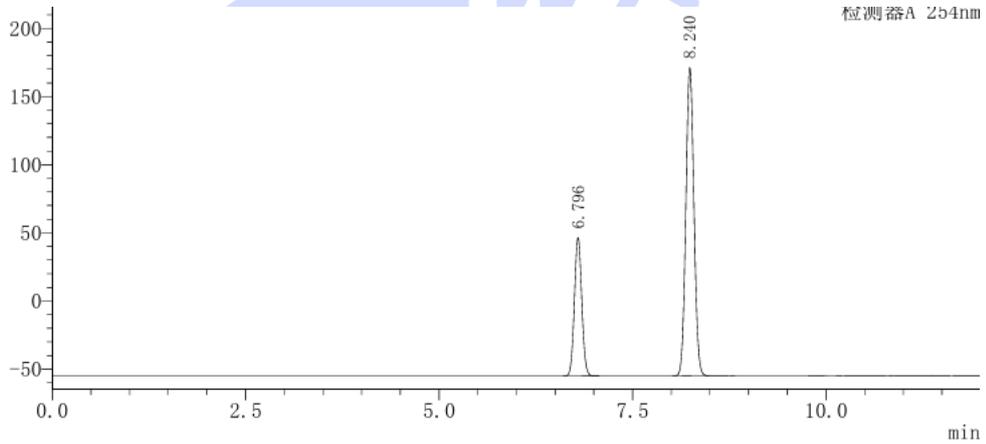## (3) Analysis spectrum of I-5 reference solution (I-5 对照溶液分析图谱):

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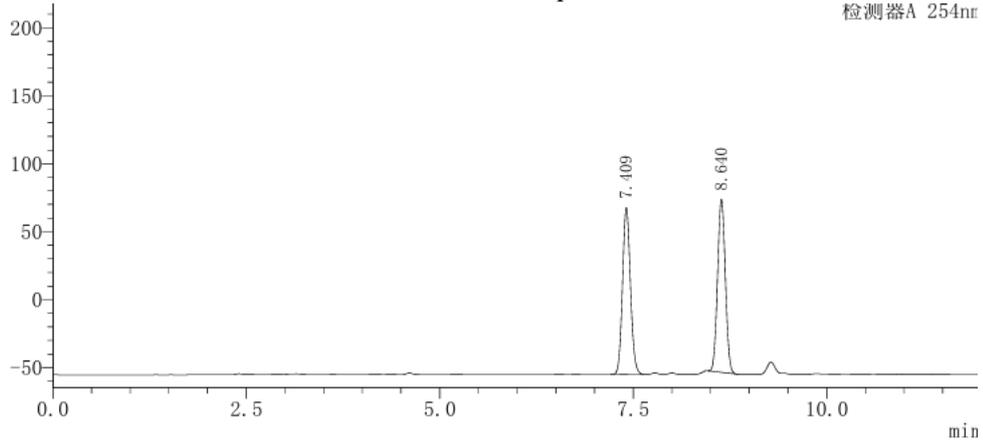
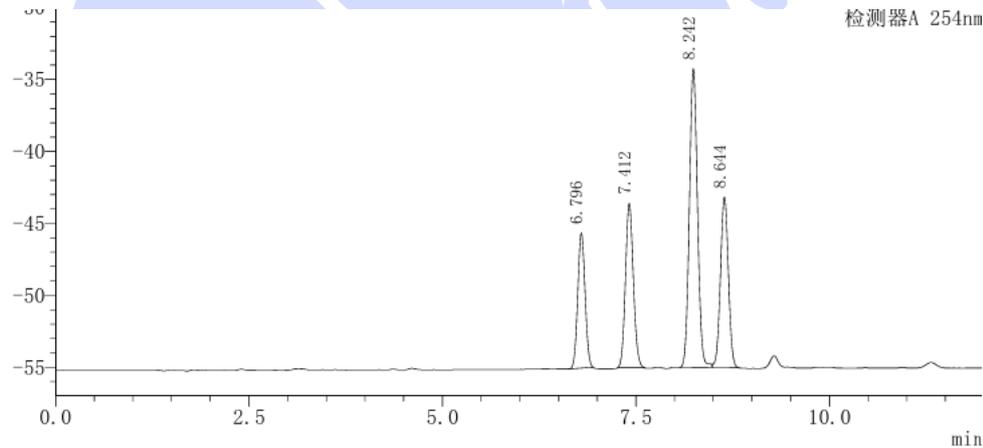
(4) Analysis spectrum of Mixed reference solution (混合对照溶液分析图谱):

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Web:www.welchmat.com

### 3. Conclusion (结论)

Using Welch Ultimate XS-C18 (4.6×150mm, 3μm) analyzed the content of Benzofurans under this chromatographic condition, all four target analytes elute within 10 minutes, and the resolution between each pair is greater than 2, that meets the customer's testing requirements.

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