Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Coccarboxylase Tetrahydrate 四水合辅羧酶	Sample properties 样品性状	Solid 固体
Date of receipt 收样日期	2025/08/14	During testing 测试期间	2025/09/02 ~ 2025/09/04
Test the ingredients and structural formula			
测试成分及结构式			
Coccarboxylase Tetrahydrate 四水合辅羧酶	Impurity Z1 杂质 Z1	Vitamin B1 维生素 B1	
Experimental Requirements			
实验要求			
Screen chromatographic columns for analyzing the system solution, with the requirement that the peak results are consistent with the typical chromatogram. 筛选色谱柱以分析系统溶液, 要求出峰结果与典型图谱一致			
Reference Method			
参考方法			
Customer's Method 客户方法			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
Water 水	Ultrapure water 超纯水	Welch 月旭	
Acetonitrile 乙腈	HPLC 色谱级	Welch 月旭	

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## 1.2. Solution preparation (溶液配制)

### 1.2.1. Preparation of mobile phase (流动相配制):

Mobile phase A: Weigh 2.40 g of Sodium dihydrogen phosphate, dissolve it in 1000 ml of pure water, adjust the pH to 6.0 with Sodium Hydroxide after full dissolution, mix well and filter to obtain the solution.

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Mobile phase A

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### 1.2.3. Test solution (供试品溶液):

Weigh 5 mg of the test sample into a 10 ml volumetric flask, dissolve it with Mobile phase A, dilute to the marked line, and mix well to obtain.

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### 1.2.4. Impurity stock solution: Take 7.5 mg of the Imp Z1 and 2.5 mg of the VB1 reference substance into a 10 ml volumetric flask, dissolve it with Mobile phase A, dilute to the marked line, and mix well to obtain.

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### 1.2.5. System suitability solution: Accurately weigh 5 mg of Cocarboxylase Tetrahydrate Reference Standard and measure 1 mL of the impurity stock solution respectively, dissolve them with a

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Tel: 400-810-6969

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Web: www.welchmat.com

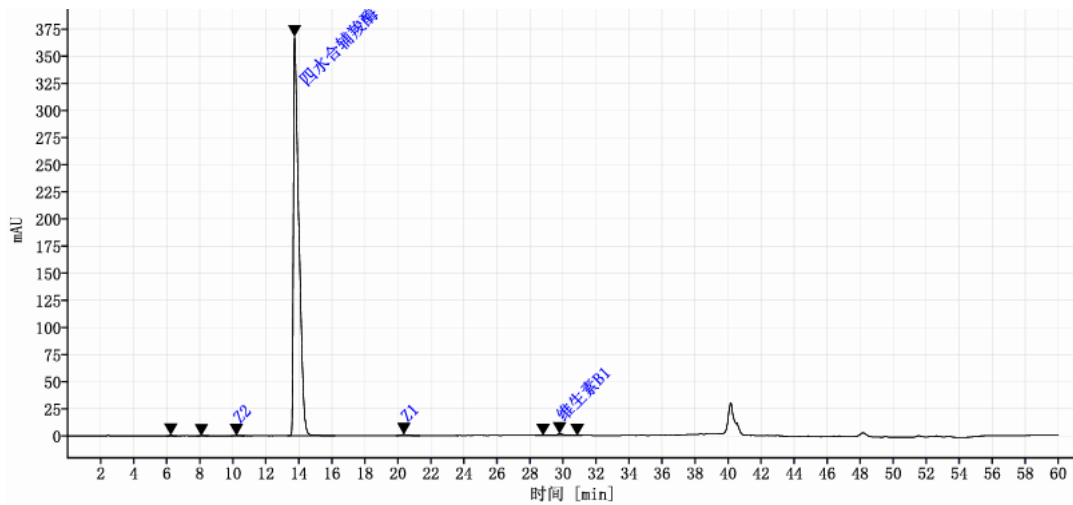
solvent, dilute to the marked volume, and mix well to obtain the solution.

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## 2. Chromatogram and data (谱图和数据)

### (1) Analysis spectrum of System suitability solution Full Range

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### (2) Analysis spectrum of System suitability solution Partial Range

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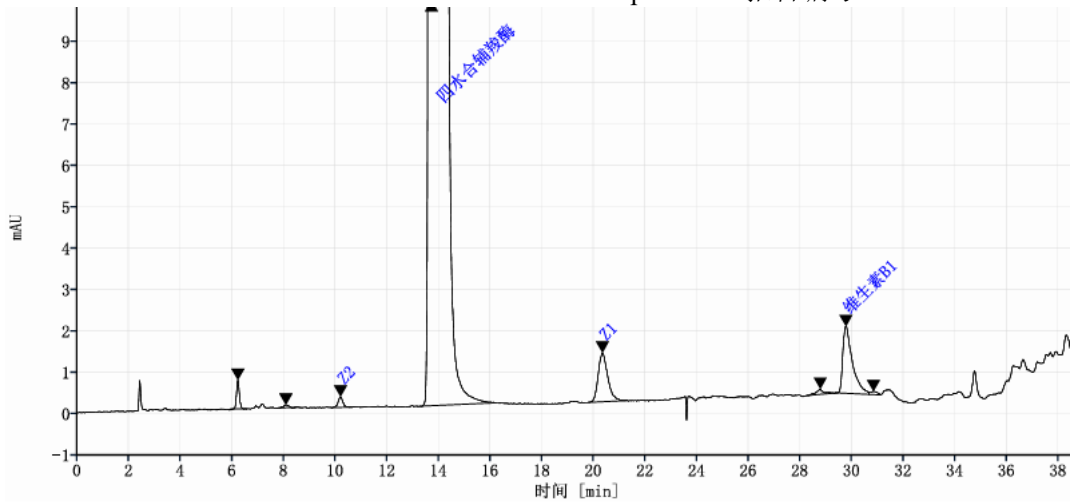
### 3. Conclusion (结论)

Using Welch Xtimate Polar-RP (4.6×250mm, 5μm) analyzed the system suitability solution under this chromatographic condition, the chromatographic results are consistent with the typical chromatogram, which meets the customer's testing requirements.

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