Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Nalfurafine Hydrochloride Related Substances 盐酸纳呋拉啡相关物质	Sample properties 样品性状	Solid 固体
Date of receipt 收样日期	2025/09/25	During testing 测试期间	2025/09/26~ 2025/09/30
Test the ingredients and structural formula			
测试成分及结构式			
<p>盐酸纳呋拉啡</p> <p>Chemical Formula: $C_{28}H_{32}N_2O_5$ Molecular Weight: 476.56</p> <p>杂质 21</p>			
Experimental Requirements			
实验要求			
<p>Develop a method to achieve effective separation of nalfurafine hydrochloride from its potential impurities (cis-impurities at positions 6a, 23, and 24; that is, Impurity 21); the requirement is that the resolution should be as high as possible.</p> <p>开发满足盐酸纳呋拉啡与其潜在杂质（6a, 23 与 24 位顺式杂质；即杂质 21）有效分离的方法；要求，分离度越大越好。</p>			

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile Phase A: Weigh 4.3 g of sodium 1-octanesulfonate, dissolve it in 1000 ml of pure water, ensure complete dissolution, adjust the pH to 6.1 with phosphoric acid, then filter and sonicate to obtain the solution.

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Methanol

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1.2.3. Test solution (测试溶液):

Test solution 1: Take an appropriate amount of nalfurafine hydrochloride, weigh it accurately, dissolve it in methanol and dilute to prepare a solution containing approximately 0.5 mg per 1 ml.

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Tel: 400-810-6969

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Web: www.welchmat.com

Test solution 2: Take an appropriate amount of Impurity 21, weigh it accurately, dissolve it in methanol and dilute to prepare a solution containing approximately 5 µg per 1 ml.

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
Test solution 3: Take appropriate amounts of nalfurafine hydrochloride and Impurity 21, weigh them accurately, dissolve them in methanol and dilute to prepare a solution containing approximately 0.5 mg of nalfurafine hydrochloride and 0.5 µg of Impurity 21 per 1 ml.

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
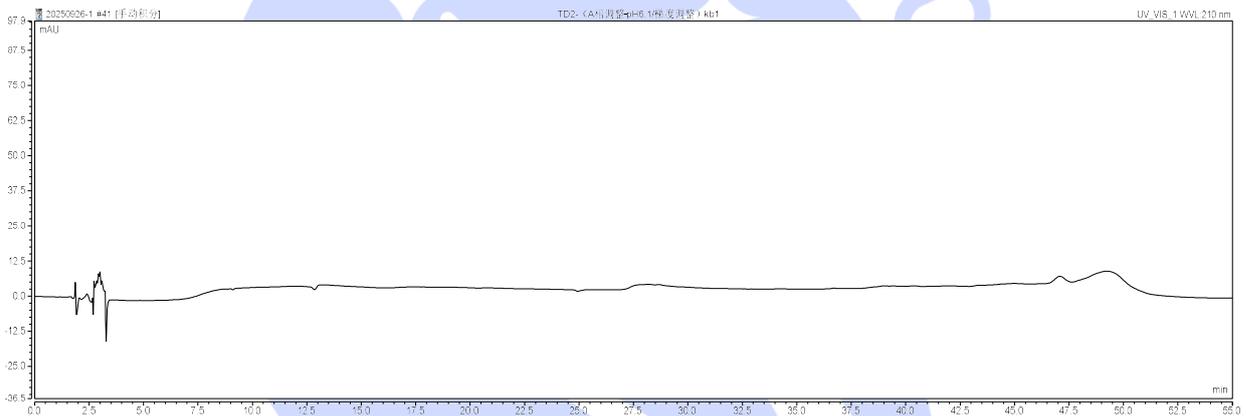
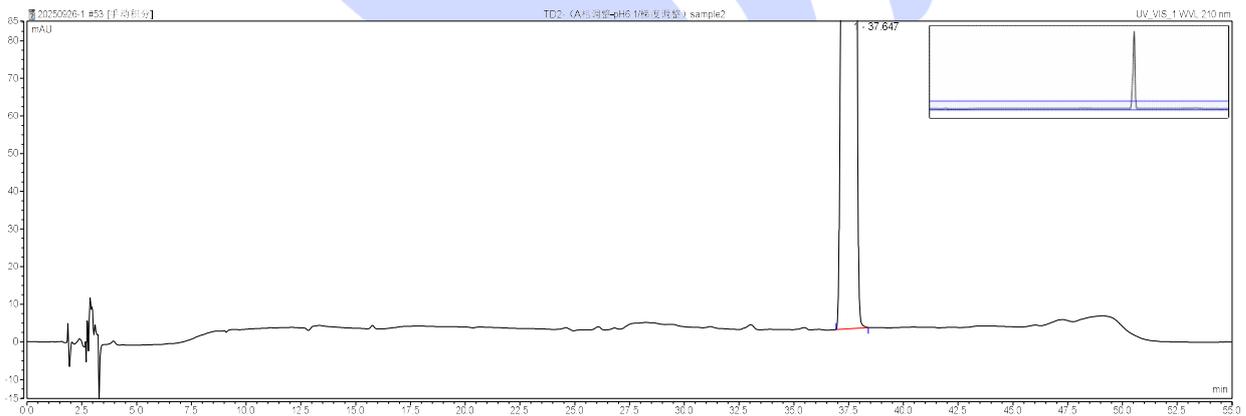
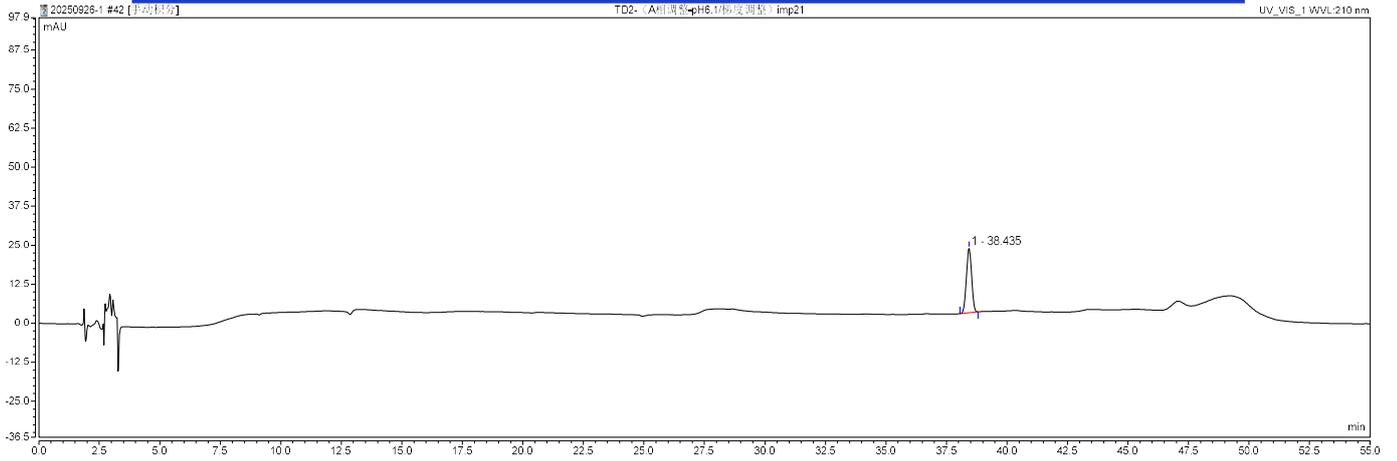
Web: www.welchmat.com

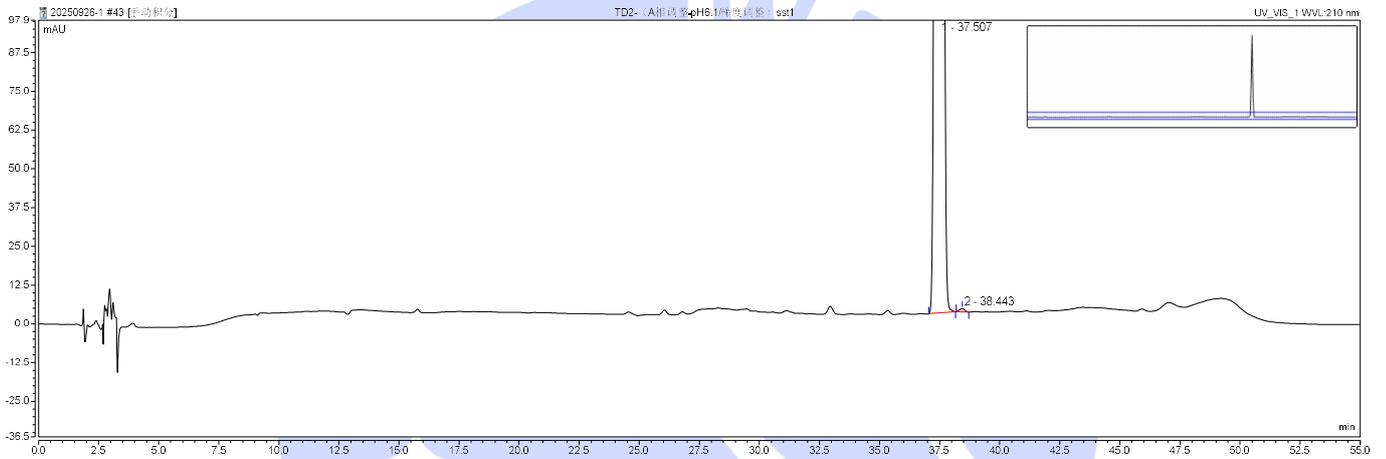(4) Analysis spectrum of Test solution 3 (测试溶液 3 分析图谱):

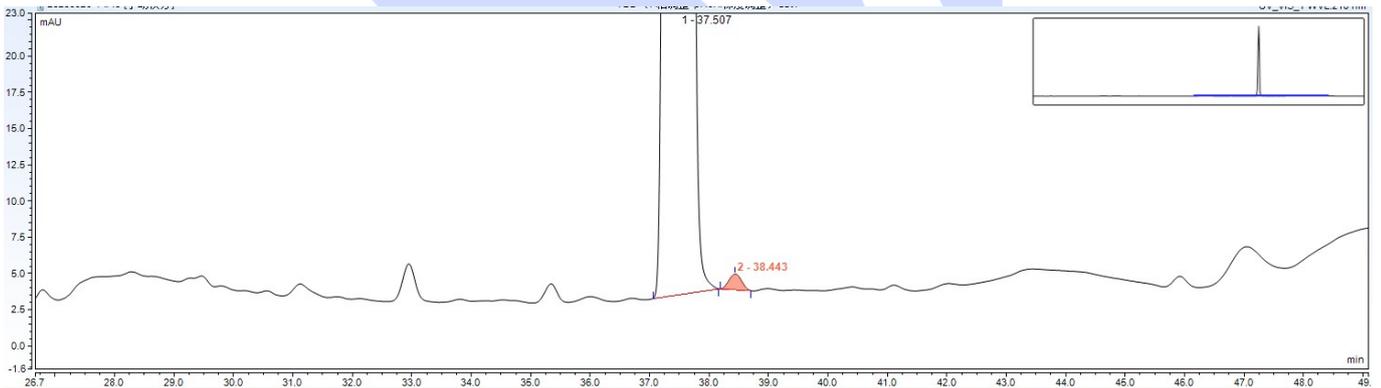(5) Chromatogram and Results of Scale-Up Analysis for Test Solution 3 (测试溶液 3 放大分析图谱及结果):

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Web: www.welchmat.com

3. Conclusion (结论)

Using the Welch Materials Xtimate-C18 chromatographic column (4.6×250 mm, 5 μm) under these chromatographic conditions, nalfurafine hydrochloride and its Impurity 21 can be effectively separated with a resolution greater than 1.5.

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