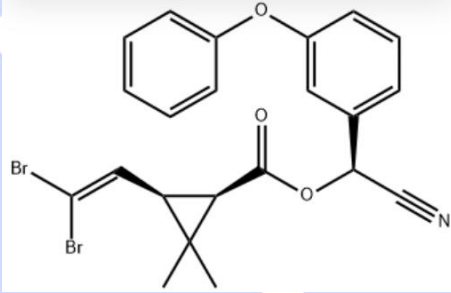
Test report

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Sample Information 样品信息			
Sample Name 样品名称	Deltamethrin and it's isomer 溴氰菊酯及异构体	Sample properties 样品性状	White solid powder 白色固体粉末
Date of receipt 收样日期	2025/07/25	During testing 测试期间	2025/08/04 ~ 08/07
Test the ingredients and structural formula 测试成分及结构式			
			
Experimental Requirements 实验要求			
<p>There is no limit on the specifications of the columns. Deltathrin and isomers are required to have a resolution of greater than 2. At present, the resolution of deltamethrin and isomers is only 1.4 under the customer method, which is relatively risky. It is best to be able to analyze four components in the same way (the mixed solution contains two components of avermectin). After deltamethrin is prepared, the longer the remaining time, the larger the isomer impurities degraded.</p> <p>柱子规格类型不限。要求溴氰菊酯和异构体的分离度大于 2。目前客户方法下溴氰菊酯和异构体分离度只有 1.4, 风险较大。最好能够同一个方法分析四个成分(混合溶液中含阿维菌素两个成分)。溴氰菊酯配制好后, 放置时间越久, 降解出的异构体杂质就越大。</p>			
Reference Method 参考方法			

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Tel:400-810-6969

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Web:www.welchmat.com

柱温	
Detector: 检测器	UV
Detection wavelength: 检测波长	245 nm
Notes 注意事项	Samples are unstable in methanol.样品在甲醇中不稳定。

1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile Phase A: Take an appropriate amount of ultrapure water and ultrasonic degassing.

Mobile Phase B: Take 800 mL of methanol, add 200 mL of methyl tert-butyl ether, mix well, and ultrasonic degassing.

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1.2.2. Test solution (供试品溶液):

Provided by customers, can be tested directly. 客户提供, 可直接测试。


1.2.3. Reference solution (对照溶液):

Weigh 0.2 mg of deltamethrin Precisely, add 1 mL of acetonitrile to dissolve.

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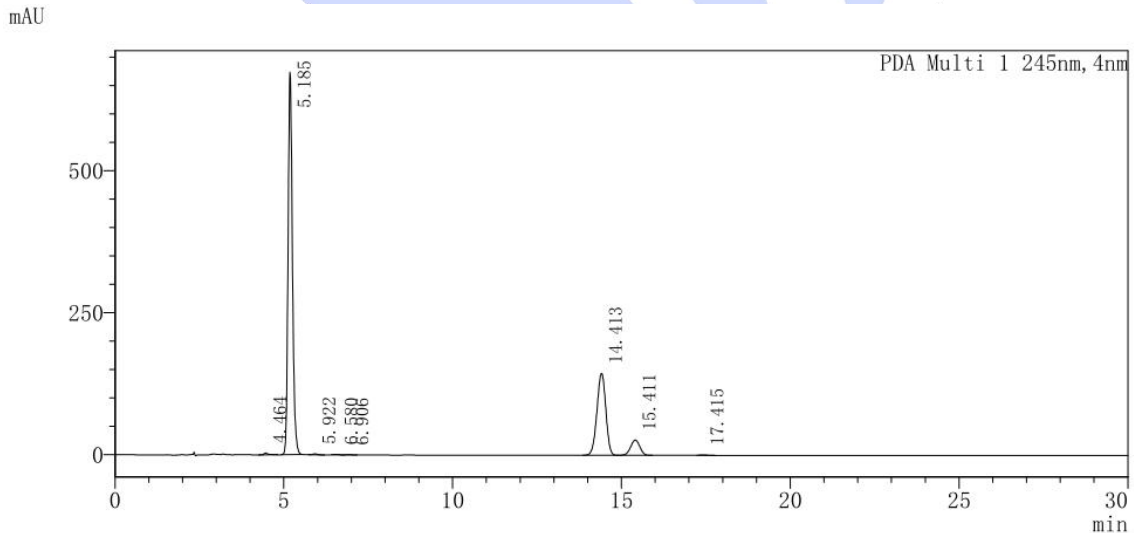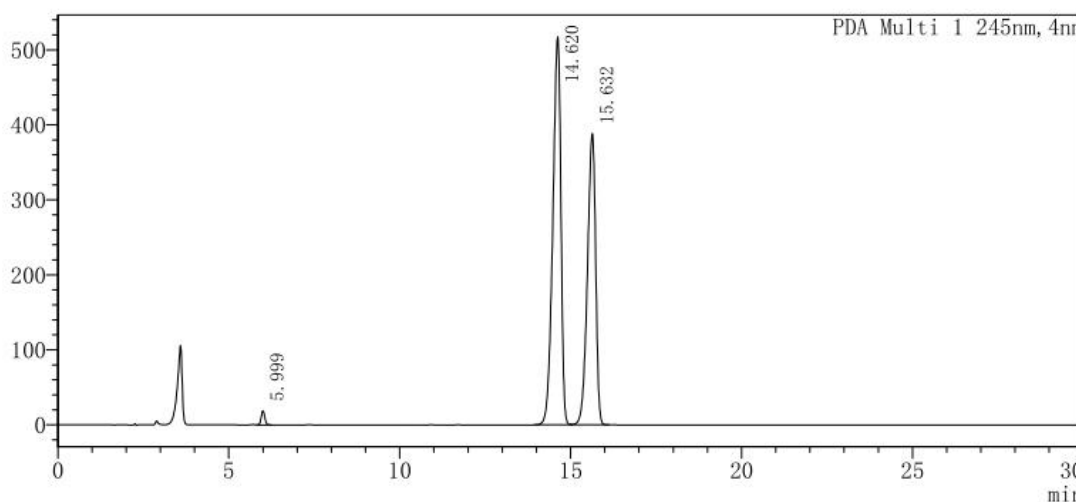
3. Conclusion (结论)

Using Welch Ultimate® XB-C18 (4.6×150mm,3µm), the detection requirements can be met under these chromatographic conditions.

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