Test report

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Sample Information 样品信息			
Sample Name 样品名称	Aminophentol oral solution 氨溴特罗口服溶液	Sample properties 样品性状	Colorless transparent liquid 无色透明液体
Date of receipt 收样日期	2025/07/25	During testing 测试期间	2025/07/25 ~ 09/08
Test the ingredients and structural formula 测试成分及结构式			
Clenbuterol 克仑特罗			
Experimental Requirements 实验要求			
The development method used UHPLC chromatography column or Boltimate chromatography column, and the peak was released within 30min 开发方法用UHPLC色谱柱或Boltimate色谱柱, 30min以内出峰			
Reference Method 参考方法			
Method provided by the client 客户方法			
Reagent information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
癸烷磺酸钠	AR	阿拉丁	
磷酸二氢钾	AR	泰坦	

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## 1.2. Solution preparation (溶液配制)

### 1.2.1. Preparation of mobile phase (流动相配制):

Mobile phase A: Weigh 3.0g sodium decanthalate and 5.0g potassium dihydrogen phosphate, add 1000ml of water to dissolve, and adjust the pH value to 3.0 with phosphoric acid.

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Mobile phase B: acetonitrile-methanol (250:150)

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### 1.2.2. Test solution (供试品溶液):

Freeze-dried preparation: add 1.1ml of ultrapure water to the test substance for dissolution, shake and mix for 1 minute, and then directly inject the sample.

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### 1.2.3. Reference solution (对照溶液):

Take an appropriate amount of clenbuterol hydrochloride reference substance, weigh it accurately, add methanol to dissolve and dilute quantitatively into a solution containing 50 $\mu$ g per 1ml, take an appropriate amount of it and add solvent to dilute quantitatively into a solution containing about 0.5 $\mu$ g per 1ml.

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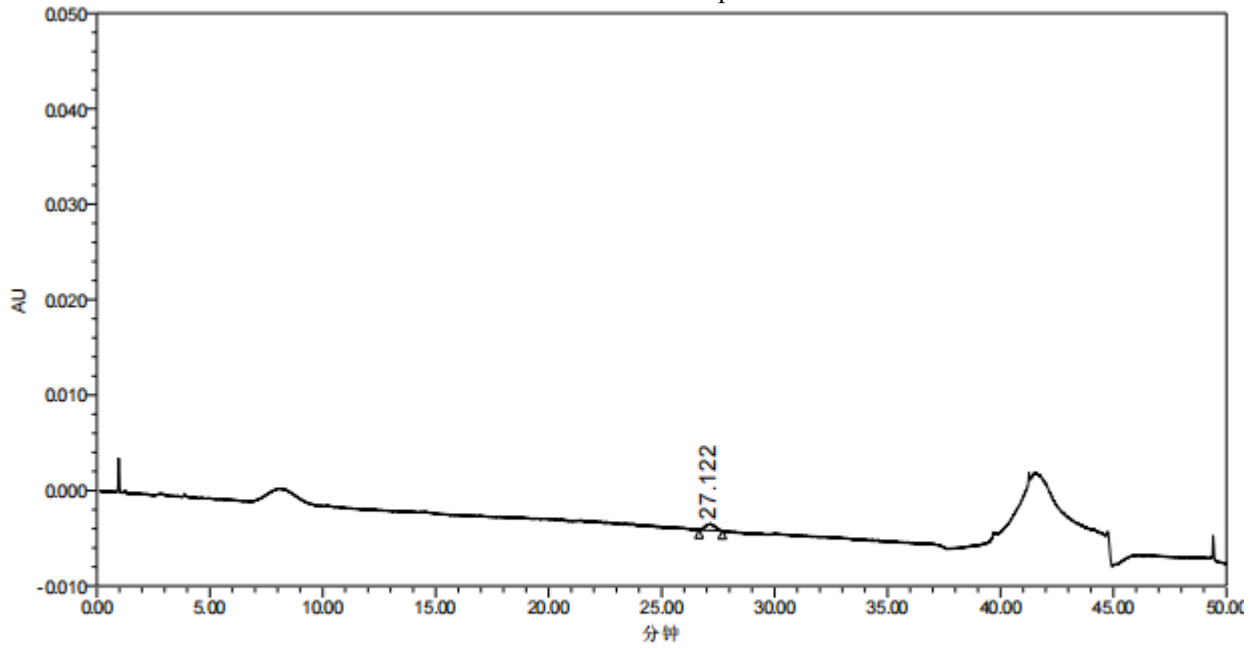
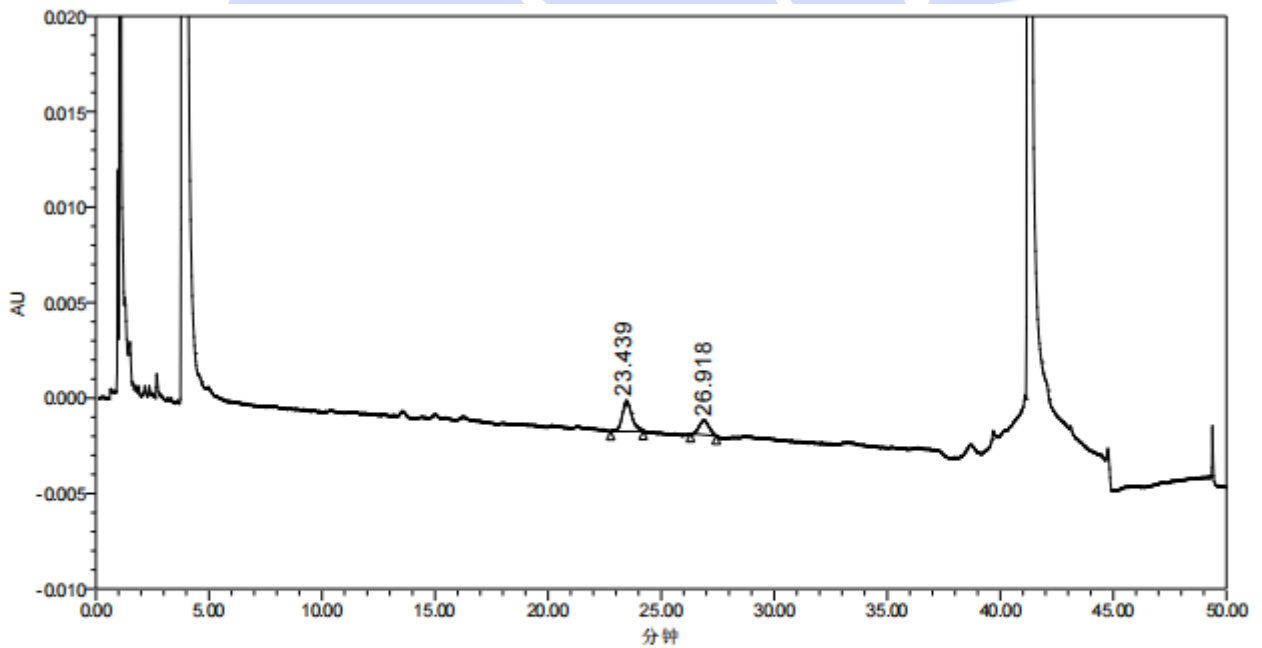
Web: www.welchmat.com

### 3. Conclusion (结论)

Using WelchUltimate<sup>®</sup> UHPLC LP-C18 (2.1×100mm,1.8μm) , Under these chromatographic conditions, the separation of reference substance and impurity peaks met the required standards. The client specified that the ambroxol concentration in the test solution should match that of the reference substance solution. During laboratory testing, no other impurities were detected in the ambroxol peak.

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