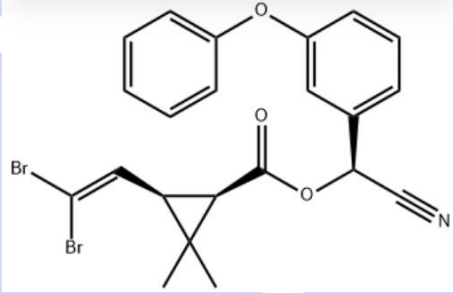
Test report

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Sample Information 样品信息			
Sample Name 样品名称	Deltamethrin and it's isomer 溴氰菊酯及异构体	Sample properties 样品性状	White solid powder 白色固体粉末
Date of receipt 收样日期	2025/07/25	During testing 测试期间	2025/09/04 ~ 09/08
Test the ingredients and structural formula 测试成分及结构式			
			
Experimental Requirements 实验要求			
<p>Appropriate methods have been developed before, but later the client hopes to redevelop a method without using methyl tert-butyl ether. Deltathrin and isomer resolution are required to be greater than 2. 此前已开发出合适方法, 但后来客户希望重新开发不使用甲基叔丁基醚的方法。要求溴氰菊酯及异构体分离度大于 2。</p>			
Reference Method 参考方法			
<p>A:Water; B:Methanol:methyl tert-butyl ether(80:20); A:B=25:75 A: 水; B: 甲醇: 甲基叔丁基醚 (80: 20); A:B=25:75</p>			
Reagent information 试剂信息			
Reagent Name	Grade	Brand	

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Web:www.welchmat.com

Report No (报告编号): 20250723-765

Mobile Phase A: Take an appropriate amount of ultrapure water and ultrasonic degassing.

Mobile Phase B1: Take 800 mL of methanol, add 200 mL of n-butanol, mix well, and ultrasonic degassing.

Mobile Phase B2: Take 800 mL of methanol, add 200 mL of Isopropanol, mix well, and ultrasonic degassing.

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1.2.2. Test solution (供试品溶液):

Provided by customers, can be tested directly. 客户提供, 可直接测试。

1.2.3. Reference solution (对照溶液):

Weigh 0.2 mg of deltamethrin Precisely, add 1 mL of acetonitrile to dissolve.

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