Test report

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| Sample Information | | | |
|---|---------------------------------|---------------------------|--------------|
| 样品信息 | | | |
| Sample Name 样品名称 | Lilarutide Injection 利拉鲁肽注射液 | Sample properties 样品性状 | Liquid 液体 |
| Date of receipt 收样日期 | 2025/08/22 | During testing 测试期间 | 2025/09/01 |
| Test the ingredients and structural formula | | | |
| 测试成分及结构式 | | | |
| Lilarutide and polymers 利拉鲁肽及聚合物 | | | |
| Experimental Requirements | | | |
| 实验要求 | | | |
| <p>The resolution in system suitability solution A shall not be less than 2.0 (calculated as the ratio of the polymer peak height to the peak-to-trough height between liraglutide and the polymer)..</p> <p>系统适用性溶液 A 中分离度应不小于 2.0 (以聚合物峰高与利拉鲁肽及聚合物之间的峰谷高之比作为分离度)</p> <p>The tailing factor of liraglutide in system suitability solution A should be between 0.8 and 2.0.</p> <p>系统适用性溶液 A 中利拉鲁肽的拖尾因子应在 0.8~2.0 之间。</p> | | | |
| Reference Method | | | |
| 参考方法 | | | |
| Customer method 客户方法 | | | |
| Reagent information | | | |
| 试剂信息 | | | |
| Reagent Name 试剂名称 | Grade 级别 | Brand 品牌 | |

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Web:www.welchmat.com

1.2.1. Preparation of mobile phase (流动相配制):

Measure 150 mL of isopropyl alcohol and 200 mL of glacial acetic acid. Add water to make up to 500 mL. Mix thoroughly, then sonicate to obtain the final solution.

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1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

Weigh approximately 1.20 g of sodium dihydrogen phosphate, dissolve in about 900 mL of water, adjust the pH to 7.5 using 1 mol/L sodium hydroxide solution, dilute to 1000 mL with water, mix thoroughly, and the solution is ready.

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1.2.3. Test solution (供试品溶液):

Accurately transfer 2.0 mL of liraglutide injection into a 25 mL volumetric flask. Add the diluent, dilute to the mark, and mix well. The solution is now ready for use.

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1.2.4. System suitability solution A (系统适用性溶液 A):

Place the liraglutide injection at 60°C for 60 hours until the polymer content reaches approximately 1.5% to 2.5%. Accurately transfer 0.84 mL of the solution into a 10 mL volumetric flask, dilute to the mark with the diluent, and mix thoroughly. The solution is now ready for use..

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1.2.5. System suitability solution B (系统适用性溶液 B):

Pipette 1.0 mL of System Suitability Solution A and 1.0 mL of Diluent, mix well, and the solution is ready..

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2. Chromatogram and data (谱图和数据)

(1) System Applicability Solution A Range Chart (系统适用性溶液 A 部分量程图谱)

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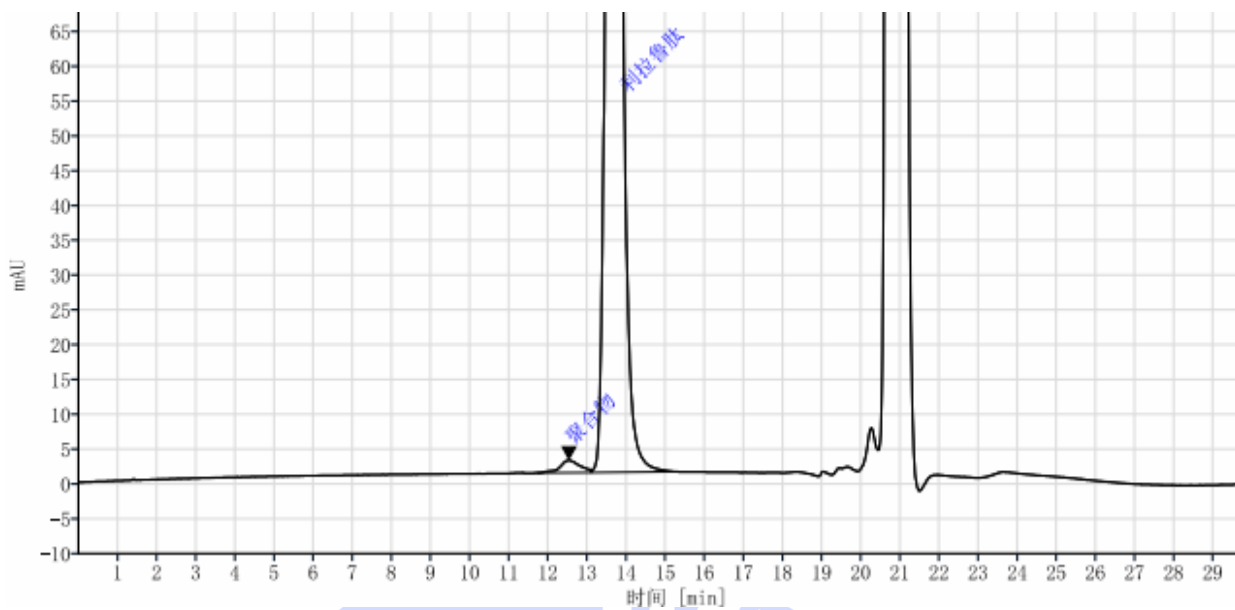
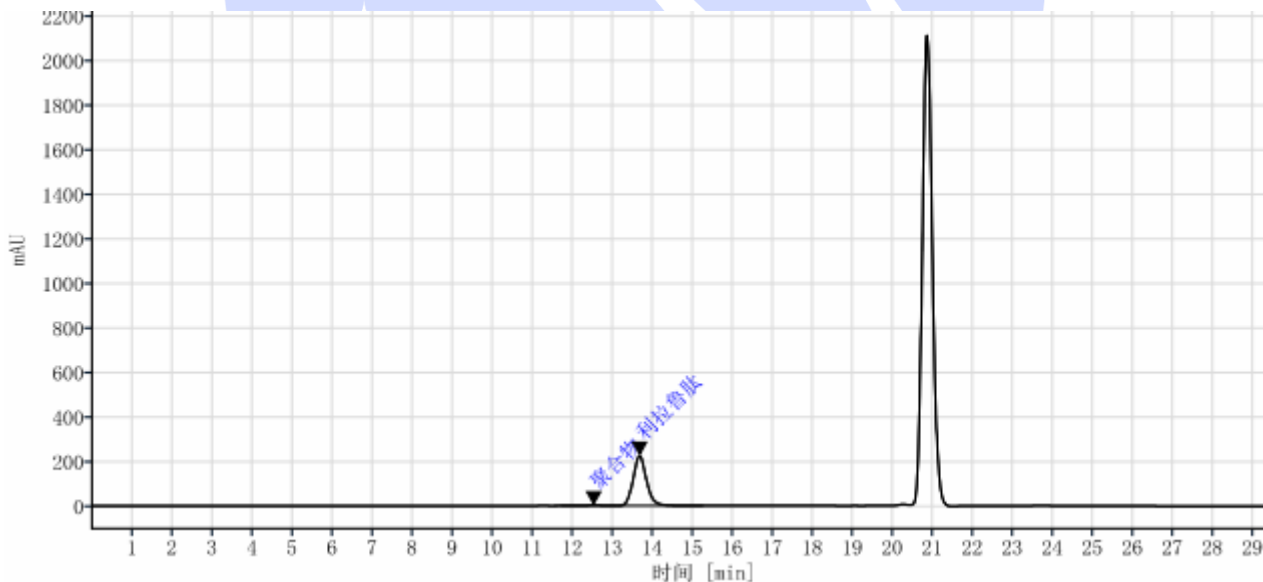
(2) System Suitability Solution A Full Range Profile (系统适用性溶液 A 全量程图谱)

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3. Conclusion (结论)

Using Welch Xtimate SEC-120A (7.8×300mm,5μm), the detection requirements can be met under these chromatographic conditions.

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