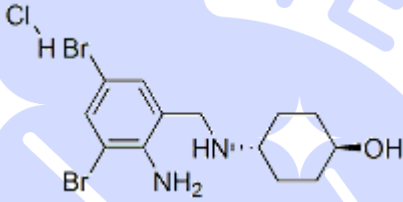
Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Ambroxol hydrochloride 盐酸氨溴索	Sample properties 样品性状	Liquid 液体
Date of receipt 收样日期	2025/11/11	During testing 测试期间	2025/11/12 ~ 2025/11/14
Test the ingredients and structural formula			
测试成分及结构式			
 <p>Ambroxol hydrochloride (盐酸氨溴索)</p>			
Experimental Requirements			
实验要求			
Analyze the test solution in accordance with the customer's method, with the requirements that the peak shape of the main component is normal and no peak splitting occurs. 按照客户方法分析供试品溶液, 要求主物质峰峰型正常, 不出现峰分叉现象			
Reference Method			
参考方法			
Customer's method 客户方法			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Article Number 货号
Water 水	Ultrapure water 超纯水	Welch 月旭	/

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile phase: Weigh 1.32 g of diammonium hydrogen phosphate and dissolve it completely in 1000 ml of purified water. Adjust the pH to 7.5 with phosphoric acid, mix well and filter. Take 480 ml of the resulting solution, mix it with 520 ml of acetonitrile, homogenize the mixture and sonicate it to obtain the final solution.

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1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

Acetonitrile-Water (1:1, v/v) (等比例乙腈水)

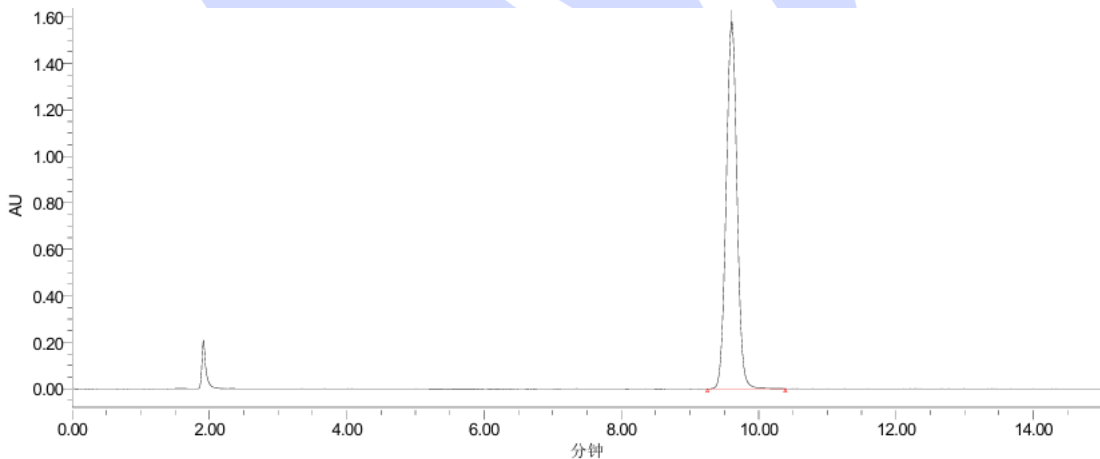
1.2.3. Test solution: Take the sample solution and dilute it 3-fold with the solvent, then mix well to obtain the final solution.

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2. Chromatogram and data (谱图和数据)

(1) Analysis spectrum of Test solution by coloum SN60250603204

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Tel: 400-810-6969

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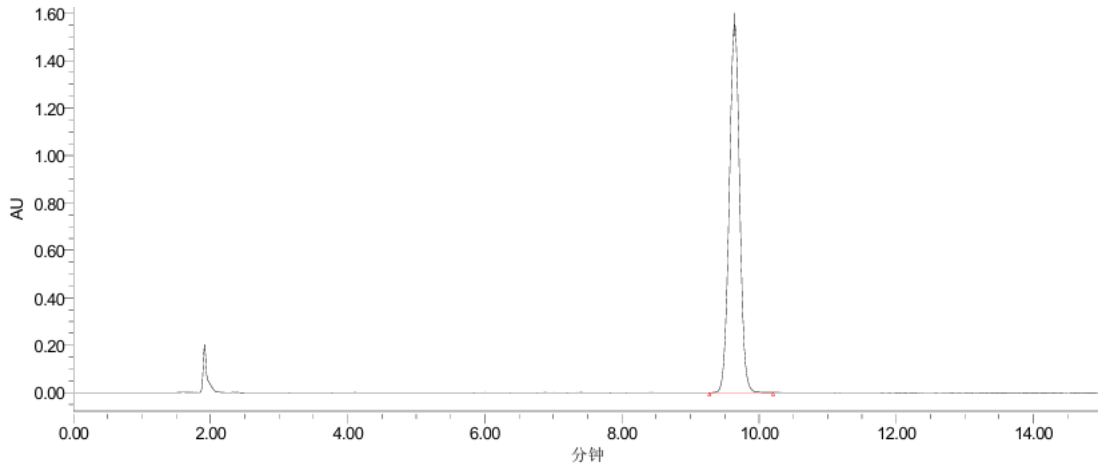
Web: www.welchmat.com

(2) Analysis spectrum of Test solution by coloum SN60250603206

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3. Conclusion (结论)

Using Welch Xtimate C18 (4.6×250mm, 5µm) analyzed test solution under this chromatographic condition, the peak of the main component eluted normally without splitting, which meets the customer's testing requirements.

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