Test report

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Sample Information 样品信息			
Sample Name 样品名称	Levofloxacin tablets 左氧氟沙星片	Sample properties 样品性状	Solid 固体
Date of receipt 收样日期	2025/11/14	During testing 测试期间	2025/11/17~11/19
Test the ingredients and structural formula 测试成分及结构式			
/			
Experimental Requirements 实验要求			
In the system suitability solution chromatogram (294nm), the resolution between the levofloxacin peak and the impurity E peak and between the levofloxacin peak and the ciprofloxacin peak should be greater than 2.0 and 2.5, respectively. 系统适用性溶液色谱中(294nm),左氧氟沙星峰与杂质E峰和左氧氟沙星峰与环丙沙星峰之间的分离度应分别大于2.0和2.5。			
Reference Method 参考方法			
Customer 客户			
Reagent information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制)

Mobile phase A: Take 4.0g of ammonium acetate and 7.0g of sodium perchlorate, add 1300mL of ultrapure water, mix well, filter, adjust pH to 2.6 with phosphoric acid solution, and mix well. Take 835mL of the above solution, add 165mL of chromatographic grade acetonitrile, shake well, and degas by ultrasonic waves.

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Mobile phase B: Take 1000mL of chromatographic grade acetonitrile, and degas by ultrasonic waves.

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1.2.2. System suitability solution (系统适用性溶液):

Provided by the customer.

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1.2.3. Test solution (供试品溶液):

Grind the text sample into powder. Take 70mg of the test ample powder, add 50mL of 0.1mol/L

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Tel: 400-810-6969

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Web: www.welchmat.com

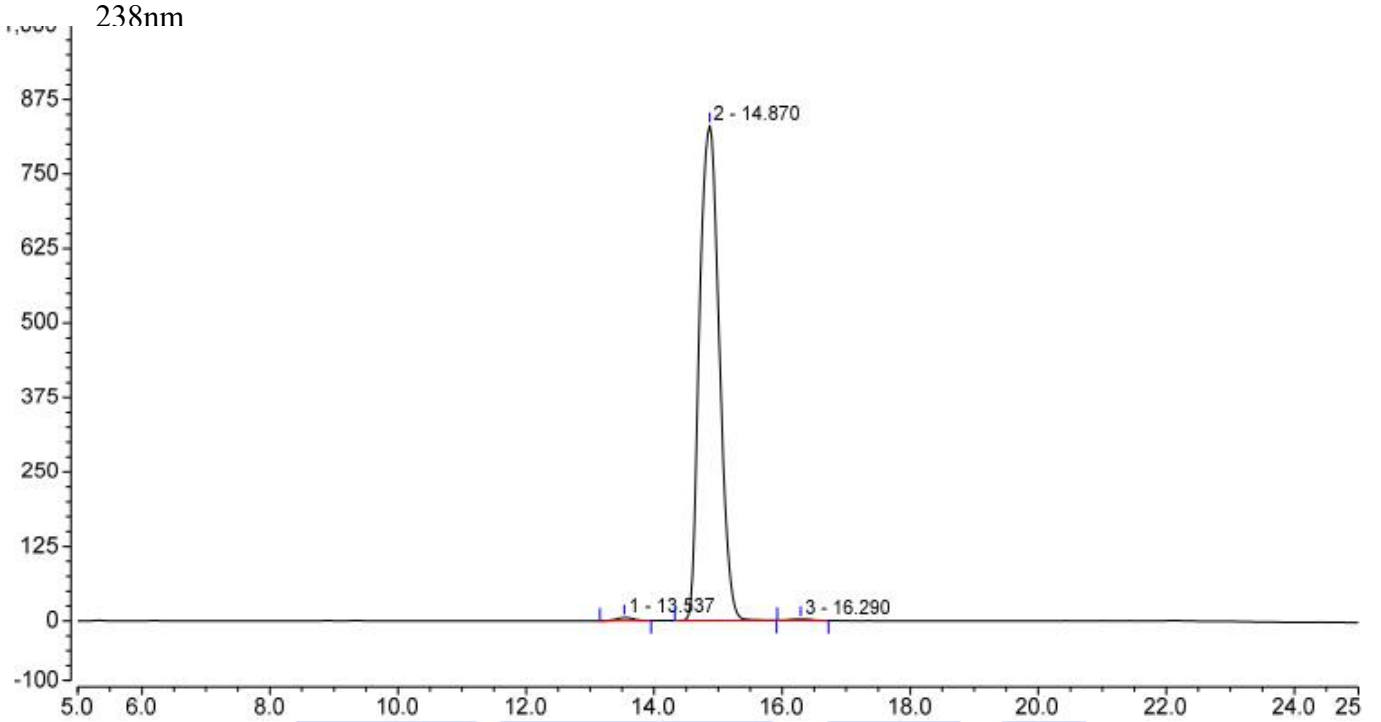
hydrochloric acid solution, mix well, and filter.

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2. Chromatogram and data (谱图和数据)

(1) System suitability solution (系统适用性溶液)

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