Test report

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Sample Information 样品信息			
Sample Name 样品名称	Glucose (15%) Electrolyte Injection 葡萄糖 (15%) 电解质注射液	Sample properties 样品性状	Colorless and transparent solution 无色透明液体
Date of receipt 收样日期	2025/11/14	During testing 测试期间	2025/12/04~12/10
Test the ingredients and structural formula 测试成分及结构式			
/			
Experimental Requirements 实验要求			
Fructose, maltose, isomaltose and maltotriose can be determined. 果糖, 麦芽糖, 异麦芽糖, 麦芽三糖能够测定。 Glucose and its impurities before and after it can be baseline separated, and blank excipients do not interfere with its determination. 葡萄糖与其前后杂质能够基线分离, 空白辅料不干扰其测定。			
Reference Method 参考方法			
No 无			
Reagent information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Item Number 货号

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Web:www.welchmat.com

Report No (报告编号): 20251107-1124-01

Take 250mL of ultrapure water, add 750mL of chromatographic grade acetonitrile, and degas by ultrasonic waves.

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1.2.2. Blank excipient solution (空白辅料溶液):

Take 200 μ L of the blank excipient solution, add 800 μ L of ultrapure water, and shake well.

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1.2.3. Mixed standard solution (混标溶液):

Take approximately 2.5mg of fructose, glucose, maltose, isomaltose, and maltotriose, respectively, dissolve each in 1mL of ultrapure water, and mix well. Take 200 μ L of each of the above five solutions, and mix well.

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1.2.4. Test solution (供试品溶液):

Take 200 μ L of the test solution, add 800 μ L of ultrapure water, and shake well.

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1.2.5. Test addition standard solution (供试加标溶液):

Take 200 μ L of the test solution, add 40 μ L of fructose solution, 100 μ L maltose solution, 400 μ L of isomaltose solution, 200 μ L of maltotriose solution, and dilute with ultrapure water to 1mL, and shake well.

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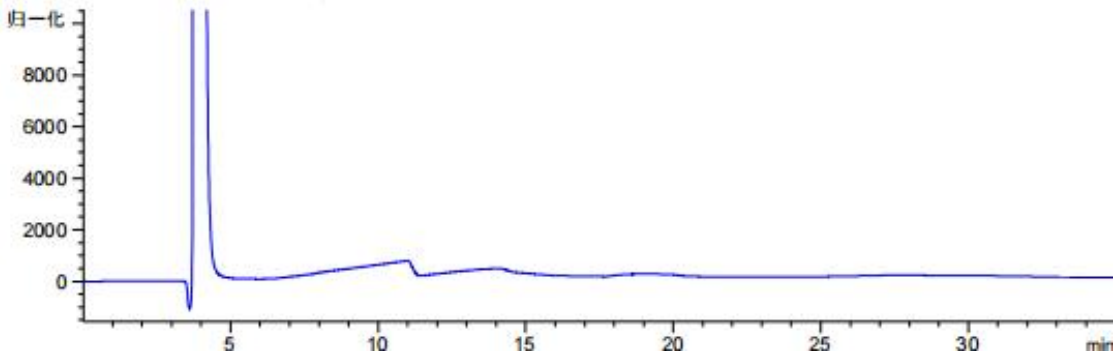
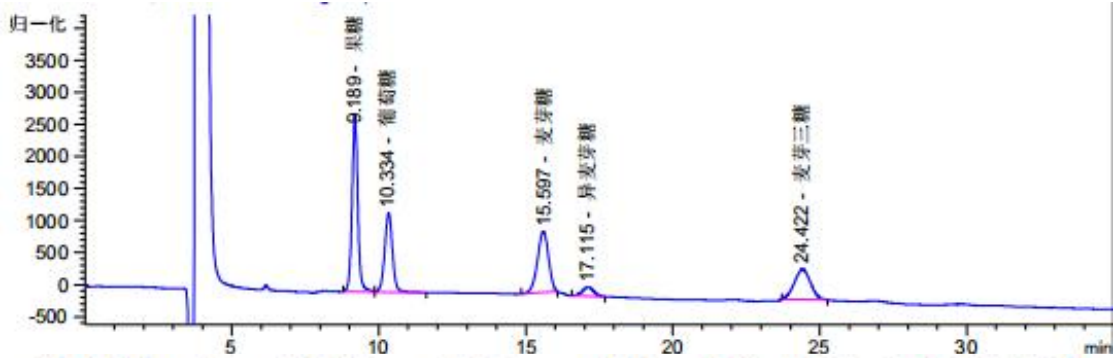
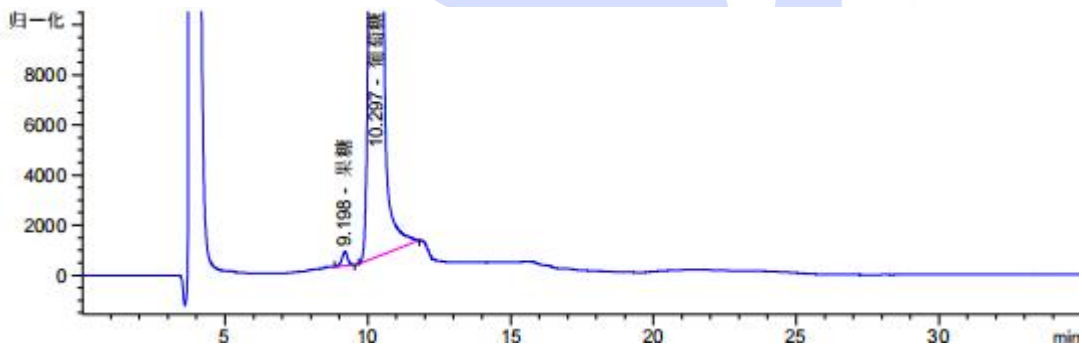
(3) Magnification view of the test solution (供试品溶液放大图)

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
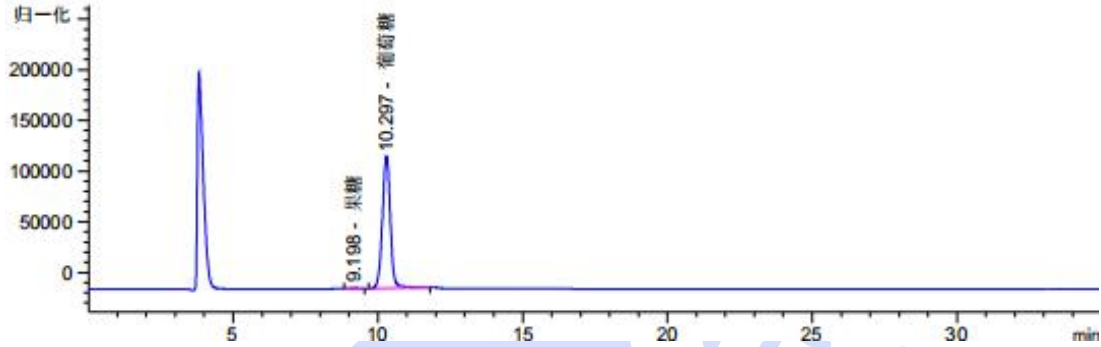
(4) Full scale of the test solution (供试品溶液满量程)

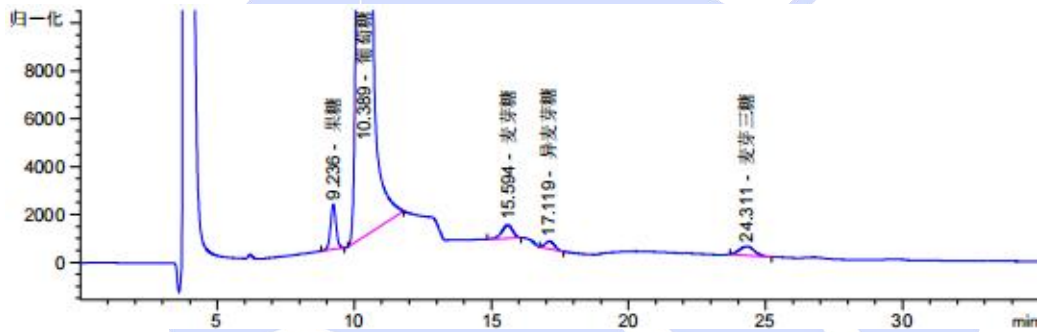
(5) Magnification view of the test addition standard solution (供试品加标溶液放大图)

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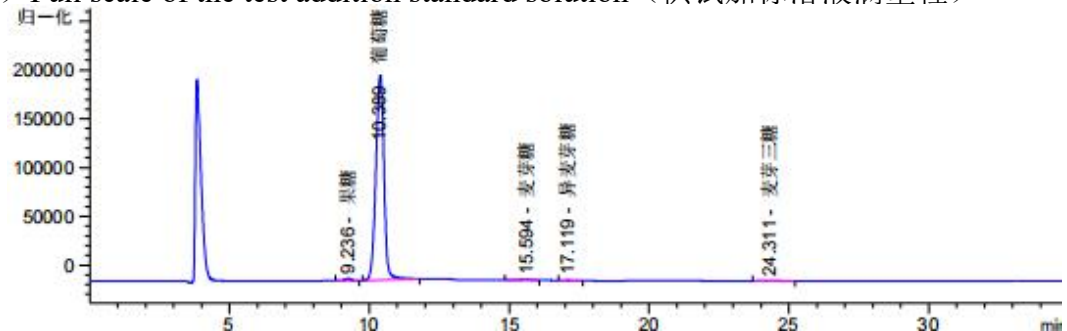(6) Full scale of the test addition standard solution (供试加标溶液满量程)

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Web:www.welchmat.com

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3. Conclusion (结论)

Using Welch Xtimate[®] NH₂ (4.6×300mm, 5μm), the detection requirements of the customer can be met under these chromatographic conditions.

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