Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Berberine Hydrochloride 盐酸小檗碱	Sample properties 样品性状	Solid 固体
Date of receipt 收样日期	2025/11/30	During testing 测试期间	2025/12/05 ~ 2025/12/08
Test the ingredients and structural formula			
测试成分及结构式			
 <p>Cl⁻</p> <p>Berberine Hydrochloride (盐酸小檗碱)</p>			
Experimental Requirements			
实验要求			
Analyze the berberine hydrochloride test solution in accordance with the customer's method, with the requirement that the main component achieves baseline separation from the adjacent subsequent impurity. 按照客户方法分析盐酸小檗碱供试品溶液, 要求主物质与后相邻杂质基线分离			
Reference Method			
参考方法			
Customer's method 客户方法			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Article Number 货号

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile phase : Weigh 1.15 g of ammonium dihydrogen phosphate and dissolve it completely in 1000 ml of purified water. Adjust the pH of the solution to 2.8 with phosphoric acid, mix well and filter. Measure 750 ml of the resulting solution, mix it with 250 ml of acetonitrile, homogenize the mixture and sonicate it to obtain the final solution.

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1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

Mobile phase (流动相)

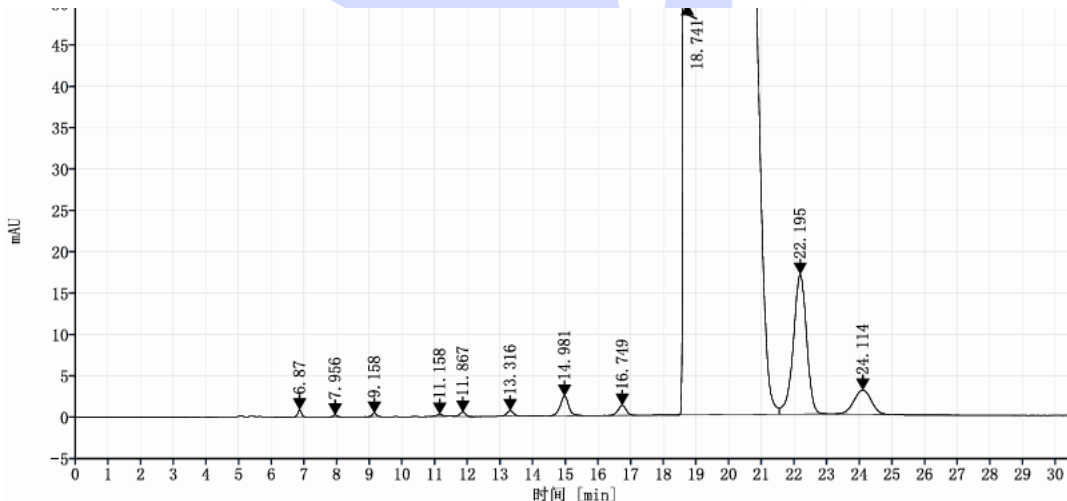
1.2.3. Test solution: Weigh 10 mg of the test substance into a 10 ml volumetric flask, dissolve and dilute it to the mark with the mobile phase, then mix well to obtain the solution.

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2. Chromatogram and data (谱图和数据)

(1) Analysis spectrum of Test solution by coloum SN60240601894

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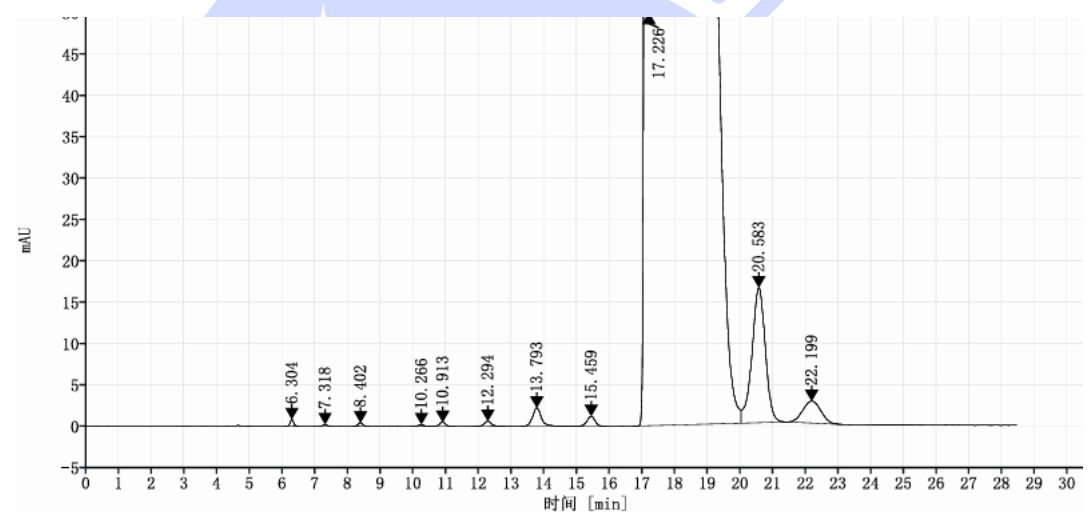
(2) Analysis spectrum of Test solution by coloum SN60250904114
(色谱柱 SN60250904114 供试品溶液分析图谱):

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Web:www.welchmat.com

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3. Conclusion (结论)

Using Welch Xtimate C18 (4.6×250mm, 5μm) analyzed test solution under this chromatographic condition, The resolution between the main component peak and the adjacent subsequent impurity peak is all greater than 1.5, which meets the customer's testing requirements.

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