Test report

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Sample Information 样品信息			
Sample Name 样品名称	Pioglitazone metformin tablets 吡格列酮二甲双胍片	Sample properties 样品性状	Solid/Liquid 固体/液体
Date of receipt 收样日期	2025/11/17	During testing 测试期间	2025/11/17~11/24
Test the ingredients and structural formula 测试成分及结构式			
/			
Experimental Requirements 实验要求			
System adaptability requirements: The resolution of metformin and p-methylacetophenone peaks shall not be less than 3, and the resolution of pioglitazone and butyl p-hydroxybenzoate peaks shall not be less than 2.5 系统适应性要求: 二甲双胍和对甲基苯乙酮峰的分离度不低于 3, 吡格列酮和对羟基苯甲酸丁酯峰的分离度不低于 2.5			
Reference Method 参考方法			
Customer 客户			
Reagent information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Item Number 货号

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制)

Mobile phase: Take 5.75g of ammonium dihydrogen phosphate, add it to 1000mL of ultrapure water, mix well, filter by suction, and degassing by ultrasonic waves; Take 500mL of the above solution and mix it with 500mL of chromatographic grade acetonitrile, shake well, then add 7.2g of sodium dodecyl sulfate, mix well, and ultrasonically degas to obtain the product.

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0.1N hydrochloric acid solution/methanol =1/1.

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1.2.3. Blank solution (空白溶液):

Take 8.5mL of hydrochloric acid, dissolve it in water and dilute it to 1000mL, and it is ready.

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Web: www.welchmat.com

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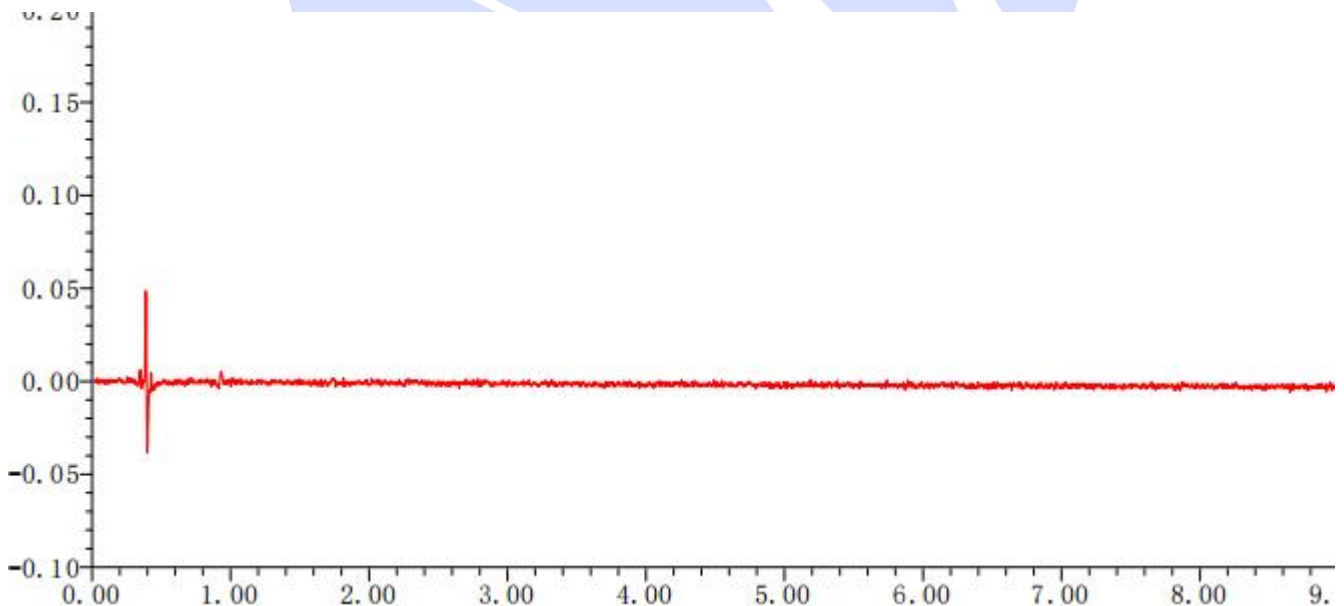
1.2.6. Test solution (供试品溶液):

Take one tablet of this product, place it in a 50mL volumetric flask, add 30mL of the diluent, ultrasonicate and shake for 60 minutes to dissolve, cool to room temperature, dilute to the mark with the diluent, and shake well. After centrifugation at 8000rpm for 5 minutes, 5 ml of the supernatant was measured and placed in a 20mL volumetric flask. The solution was diluted to the mark with the diluent and shaken well. Measure 5mL and place it in a 50mL volumetric flask. Dilute it to the mark with 0.1N hydrochloric acid solution and shake well. After filtration through a 0.45 μ m filter membrane, the filtrate is taken and the result is obtained.

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Tel: 400-810-6969

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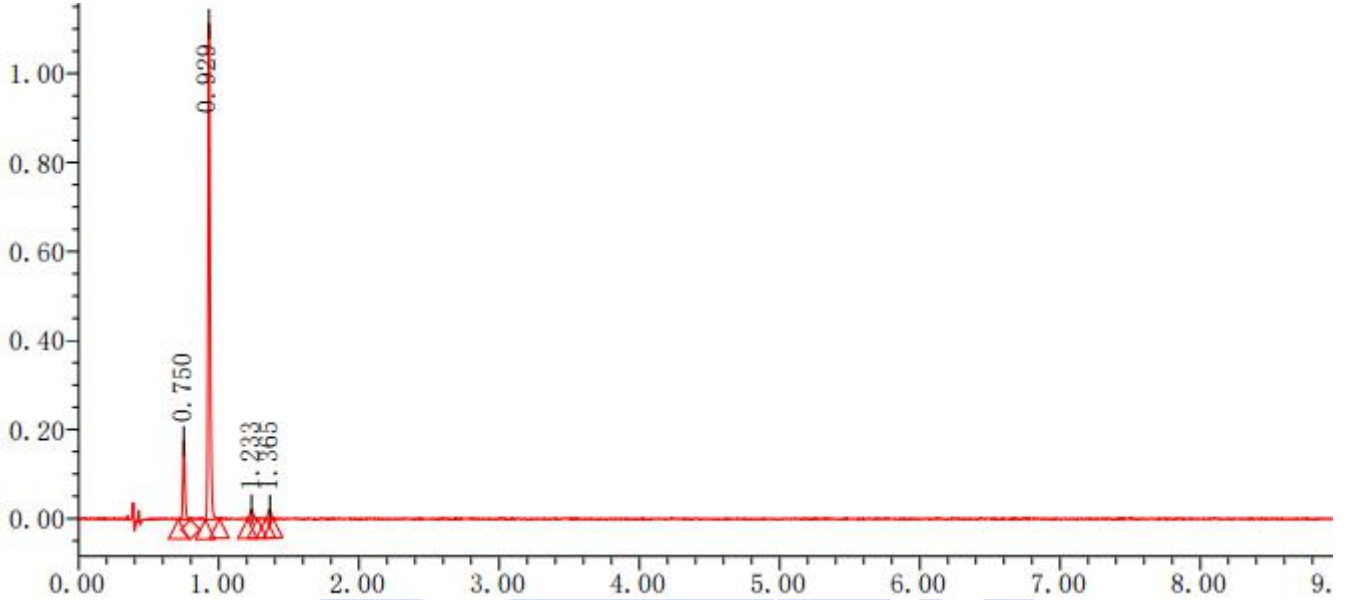
Web: www.welchmat.com


(2) System adaptive solution diagram (系统适应性图)

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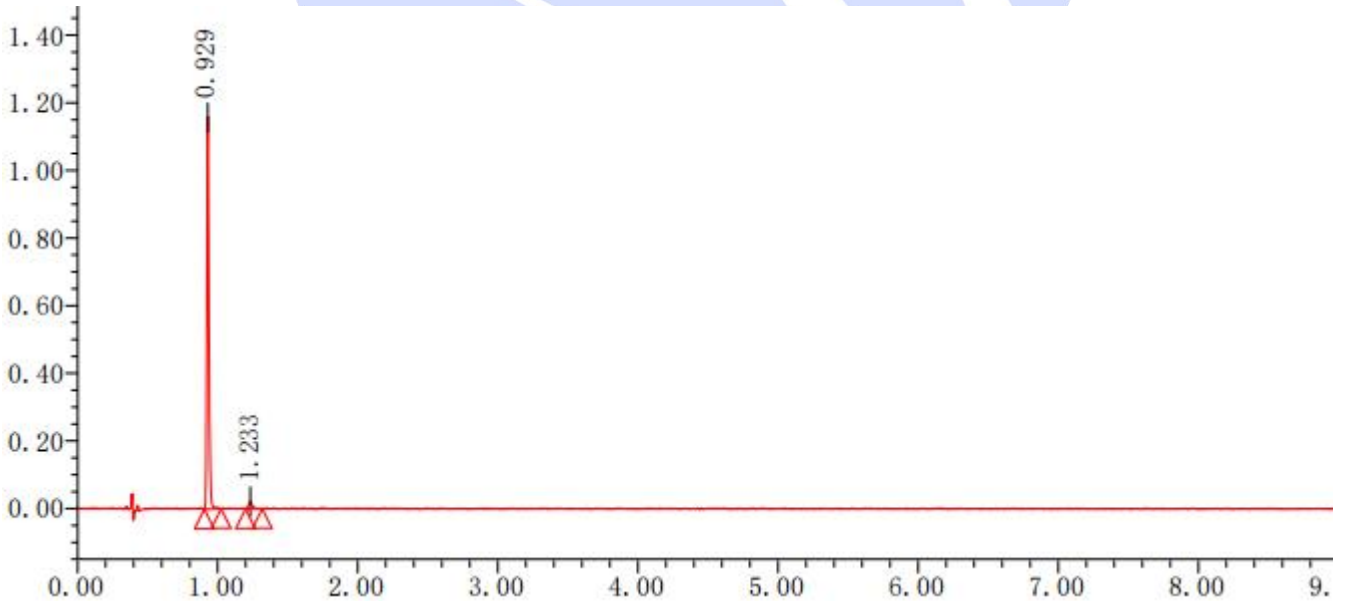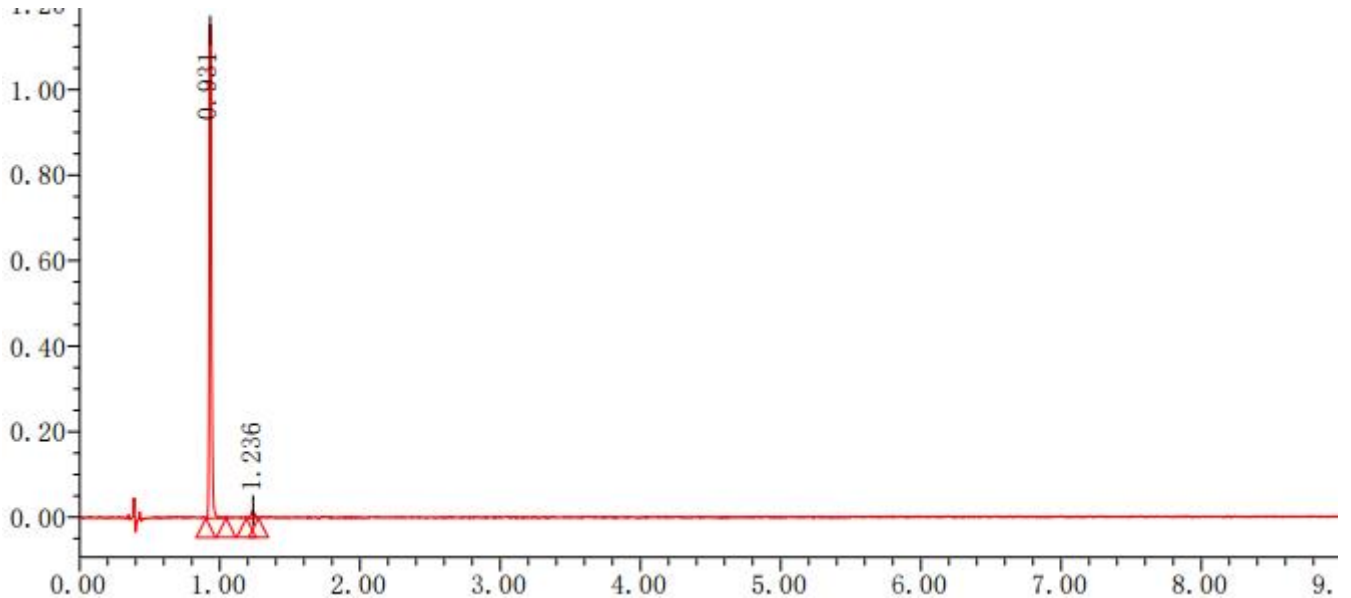
(4) Diagram of the test solution (供试品溶液图)

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3. Conclusion (结论)

Using Welch Xtimate® UHPLC C18 (2.1×100mm, 1.8µm), the detection requirements of the customer can be met under these chromatographic conditions.

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