Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Iopromide and its Chloro Impurities 碘普罗胺及其氯代杂质	Sample properties 样品性状	Liquid 液体
Date of receipt 收样日期	2025/11/01	During testing 测试期间	2025/11/05 ~ 2025/11/10
Test the ingredients and structural formula			
测试成分及结构式			
			
Iopromide (碘普罗胺)	Chloro Impuritie 1 (氯代杂质 1)	Chloro Impuritie 2 (氯代杂质 2)	
Experimental Requirements			
实验要求			
Develop a method to analyze Iopromide and its Chloro Impurities, requiring the separation between Iopromide and the impurities to meet the customer's requirements. 开发方法分析碘普罗胺及其氯代杂质，要求碘普罗胺与杂质的分离满足客户要求			
Reference Method			
参考方法			
Customer's method 客户方法			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Article Number 货号
Water 水	Ultrapure water 超纯水	Welch 月旭	/

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile Phase: Take 0.68g of Potassium Dihydrogen Phosphate into 1000ml of pure water. After thorough dissolution, adjust the pH to 6.8 with a Sodium Hydroxide solution. After mixing, take 950ml of the resulting solution and mix it with 50ml of methanol. Mix well and sonicate to obtain the final solution.

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1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

Provided by the customer (客户提供)

1.2.3. Iopromide, Impurity 1, and Impurity 2 reference solution: Provided by the customer, ready for direct analysis.

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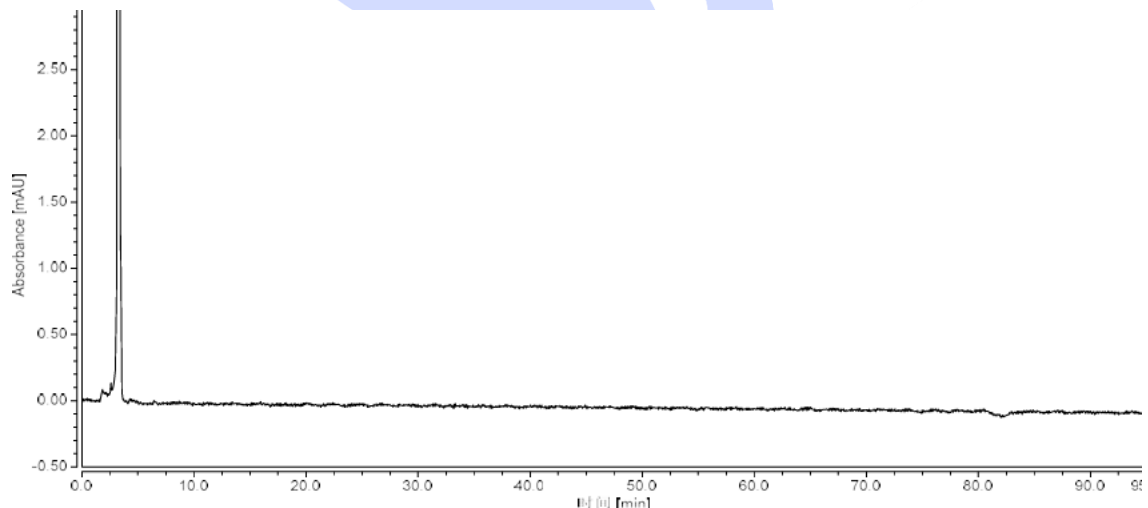
1.2.4. Mixed reference solution: Prepare by mixing each reference solution in equal proportions.

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Tel: 400-810-6969

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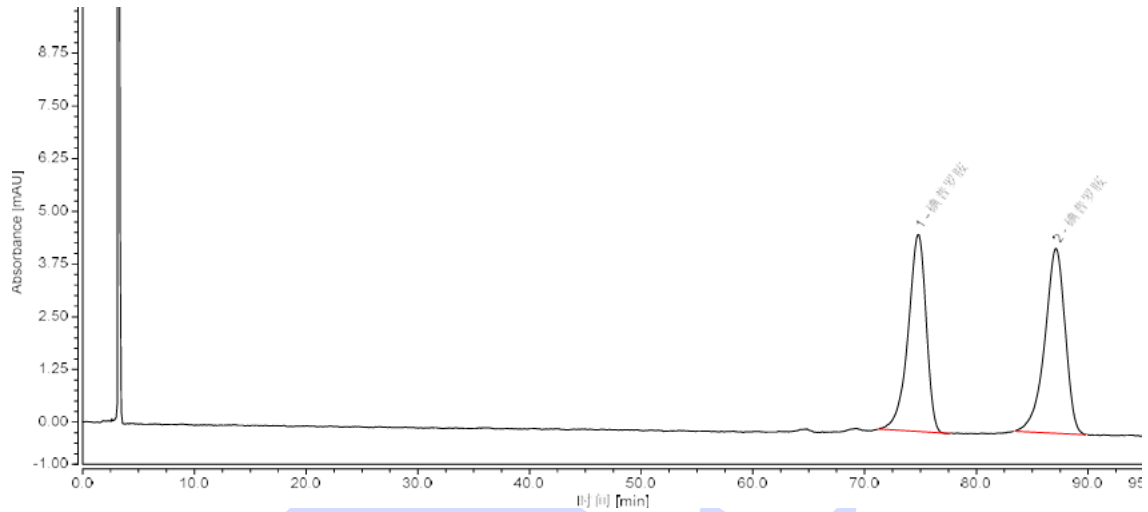
Web: www.welchmat.com


(2) Analysis spectrum of API reference solution (碘普罗胺对照溶液分析图谱):

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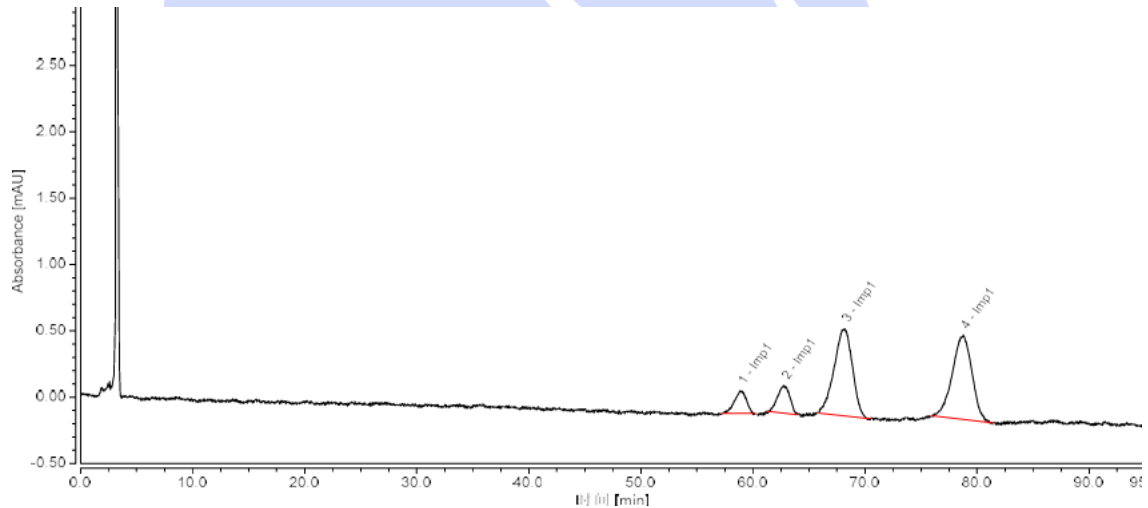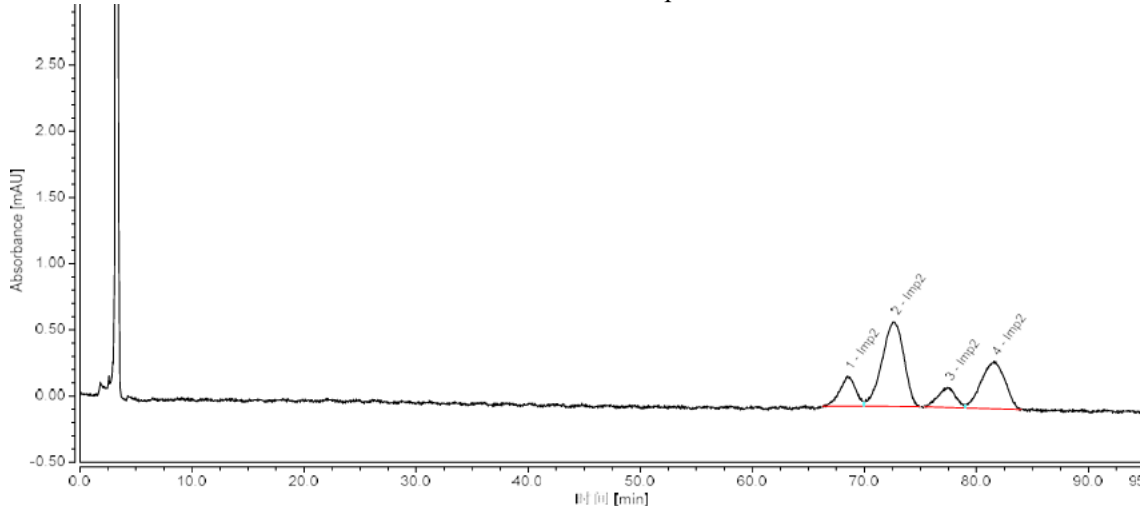
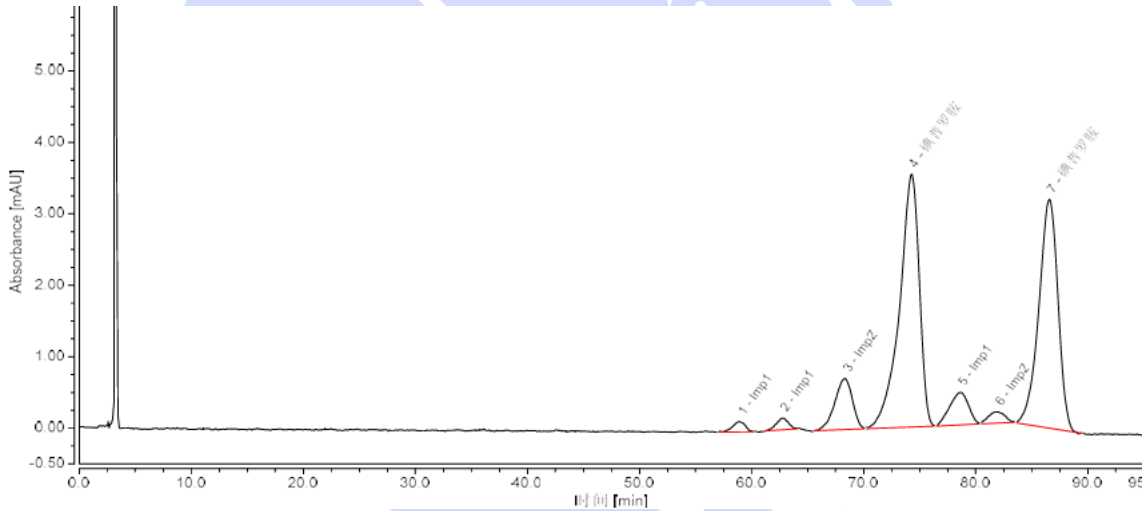
(5) Analysis spectrum of Mixed reference solution (混合对照溶液分析图谱):

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Web:www.welchmat.com

3. Conclusion (结论)

Using Welch Xtimate Phenyl-Hexyl (4.6×250mm, 5μm) analyzed Iopromide and its Chloro Impurities under this chromatographic condition , the results meets the customer's detection requirements

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