Test report

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Sample Information 样品信息			
Sample Name 样品名称	Fluvastatin Sodium Capsules 氟伐他汀钠胶囊	Sample properties 样品性状	White powder 白色粉末
Date of receipt 收样日期	2026/01/14	During testing 测试期间	2026/01/15~01/26
Test the ingredients and structural formula 测试成分及结构式			
/			
Experimental Requirements 实验要求			
<p>Screen the chromatographic column. The retention time of the fluvastatin peak in the chromatogram of the system suitability solution should be approximately 10 minutes, and the resolution between the fluvastatin peak and the impurity I peak should be greater than 1.4.</p> <p>筛选色谱柱, 要求系统适用性溶液色谱图中氟伐他汀峰保留时间约为 10 分钟, 氟伐他汀峰与杂质 I 峰之间的分离度应大于 1.4。</p> <p>In the chromatogram of the sensitive solution, the signal-to-noise ratio of the peak height of fluvastatin should be no less than 10.</p> <p>灵敏度溶液色谱图中氟伐他汀峰峰高的信噪比应不小于 10。</p>			
Reference Method 参考方法			
Customer 客户			
Reagent information			

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## 1.2. Solution preparation (溶液配制)

### 1.2.1. Preparation of mobile phase (流动相配制)

**Solution I:** Take 40 mL of 25% tetramethylammonium hydroxide solution, dilute to 1000 mL with ultrapure water, mix well, adjust the pH to 7.2 with phosphoric acid solution, and filter under suction to obtain the desired solution.

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**Solution II:** Take 400mL of chromatographic grade acetonitrile, add it to 800mL of

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Tel: 400-810-6969

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Web: www.welchmat.com

chromatographic grade methanol, and mix well.

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Mobile phase A: Take 125mL of solution II, add it to 800mL of solution I, mix well, and degas by ultrasonic waves.

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Mobile phase B: Take 125mL of solution I, add it to 800mL of solution II, mix well, and degas by ultrasonic waves.

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Diluent: Take 46mL of solution II, add it to 54mL of solution I, and mix well.

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## 1.2.2. System suitability solution (系统适用性溶液):

Take 4mg of flurvastatin sodium, add 2mL of diluent, mix well, add 200 $\mu$ L of 1mol/L hydrochloric acid solution, let stand for 30 minutes, add 100 $\mu$ L of 2mol/L sodium hydroxide solution, mix well, add diluent to 10mL, and shake well.

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## 1.2.3. Test solution (供试品溶液):

Take 45mg of the test sample powder, add 10mL of diluent, mix well, and filter.

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## 1.2.4. Sensitive solution (灵敏度溶液):

Take 5 $\mu$ L of the test solution, add diluent to 1mL, mix well. Take the above solution, add 900 $\mu$ L of diluent, and mix well.

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

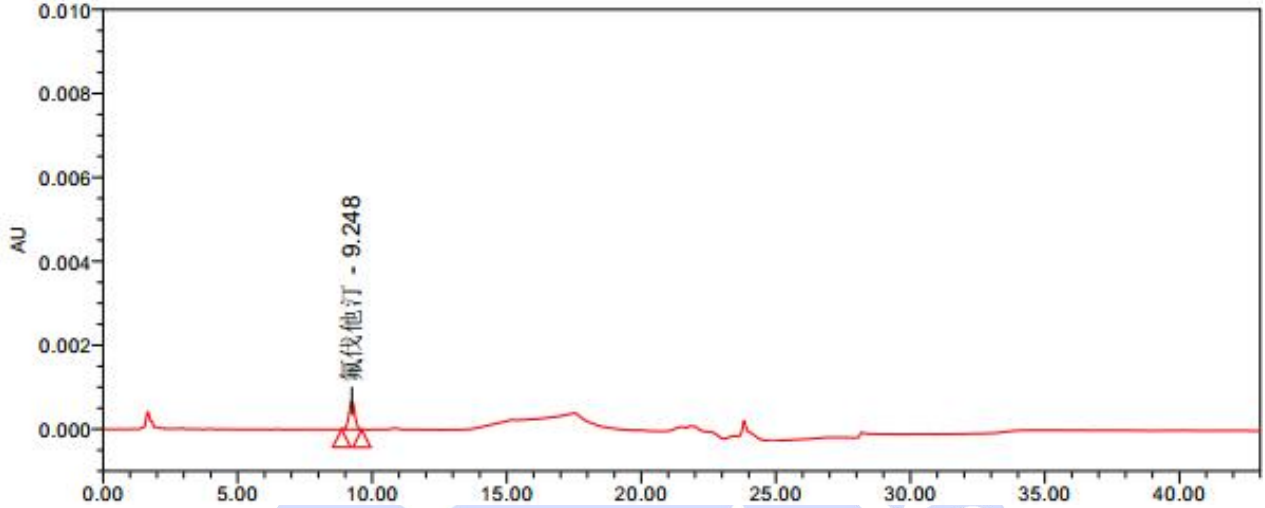
### (2) System suitability solution (系统适用性溶液图)

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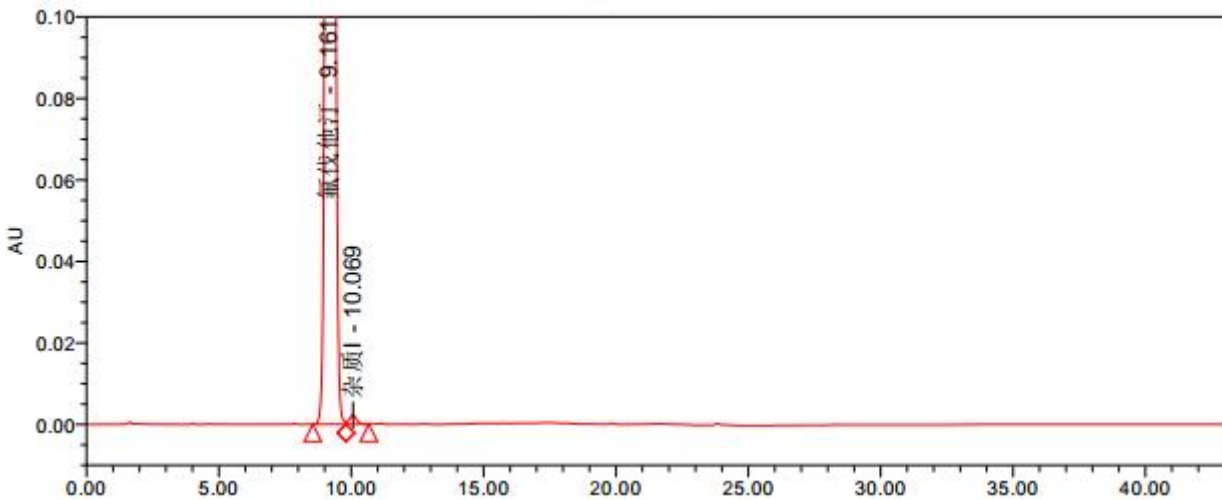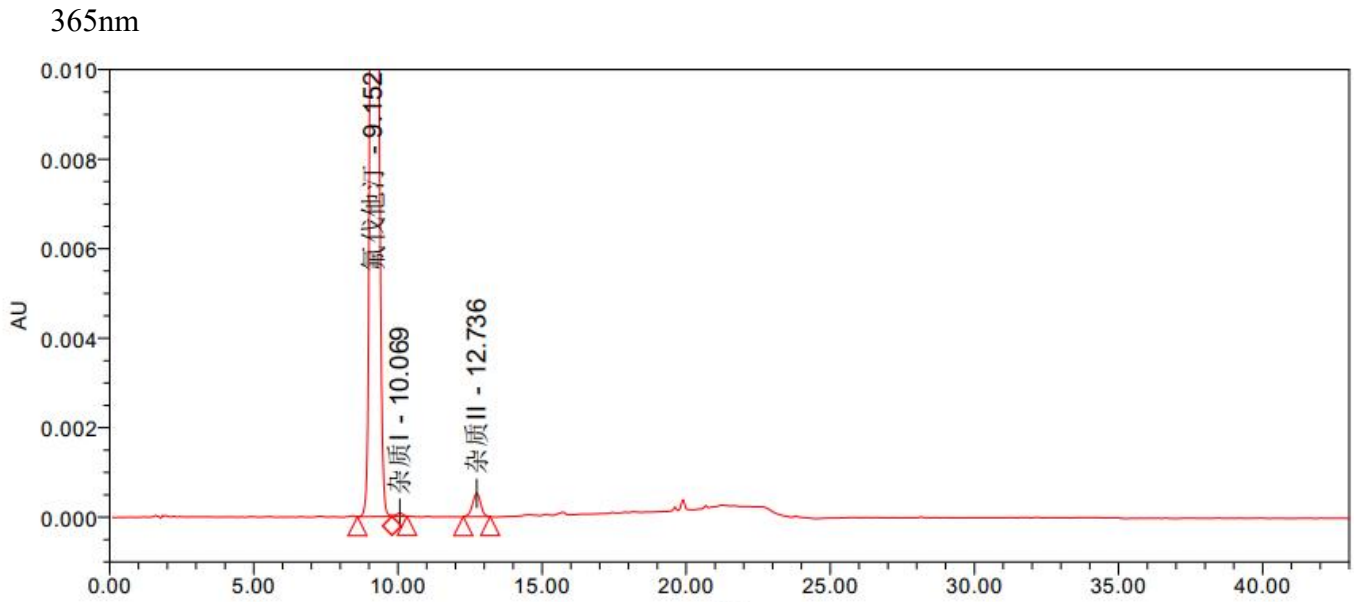
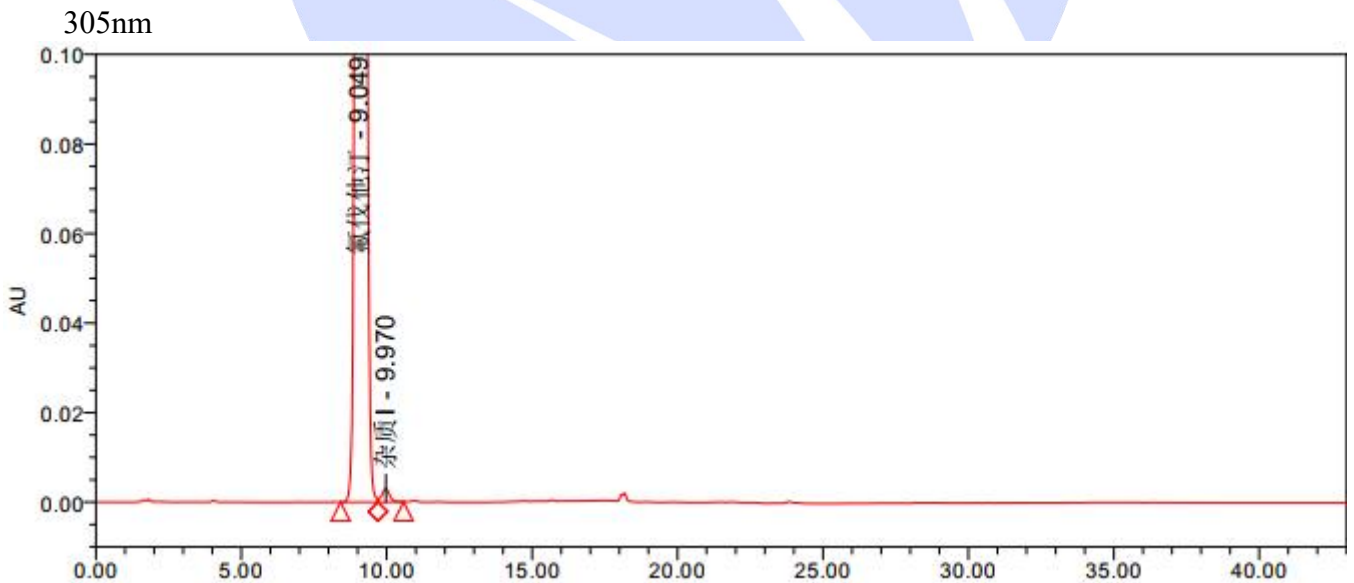
(3) Magnification view of the test solution (供试品溶液放大图)

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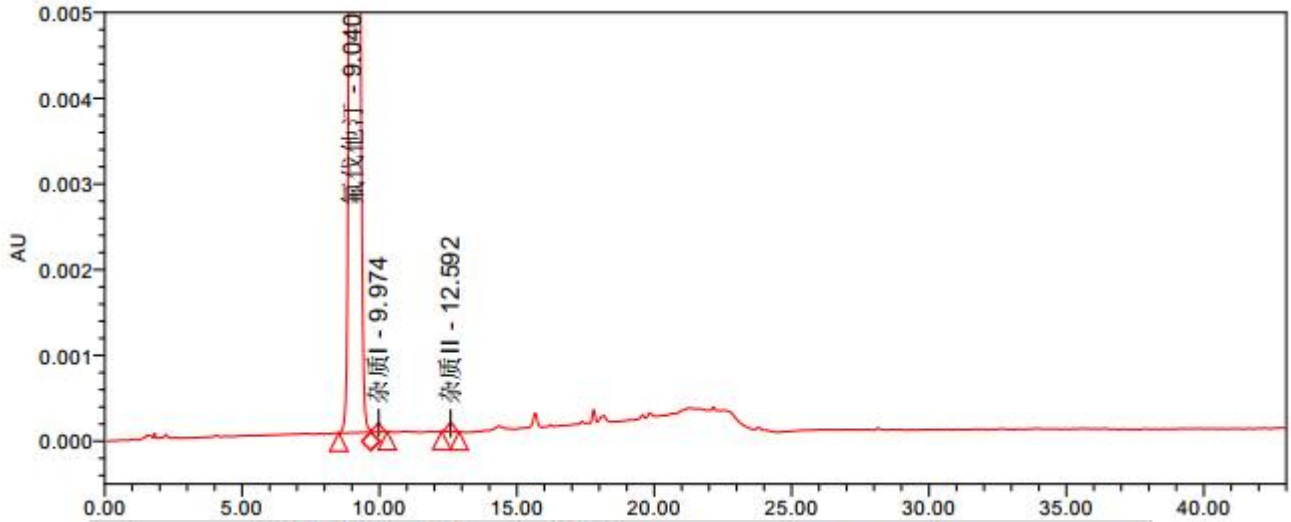
(4) Full-scale diagram of the test solution (供试品溶液满量程图)

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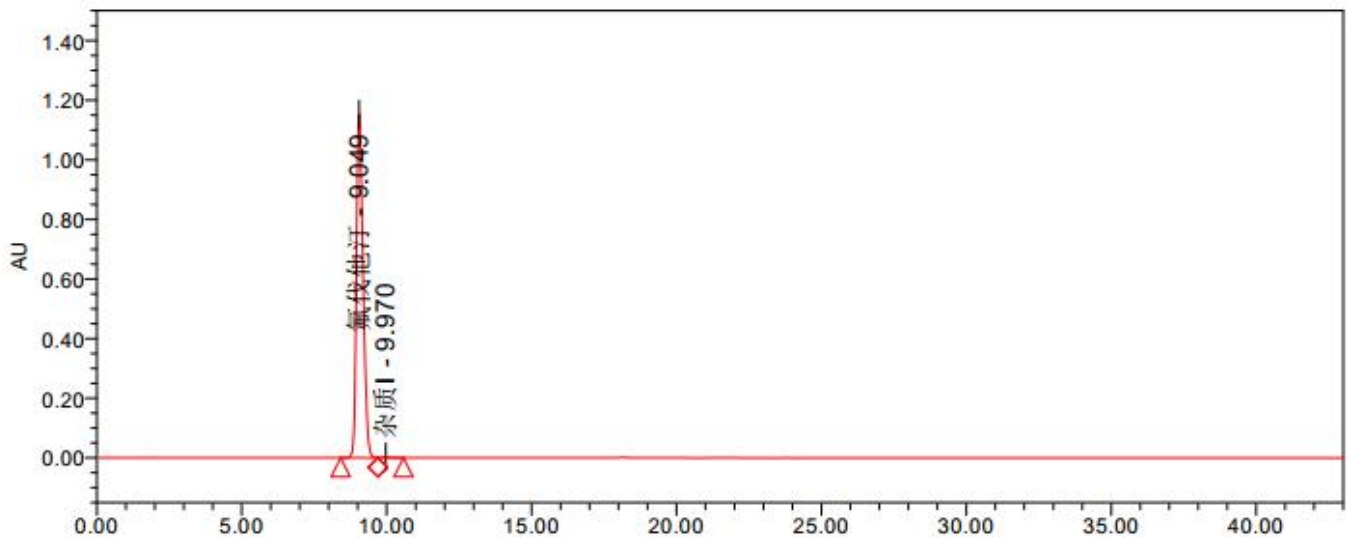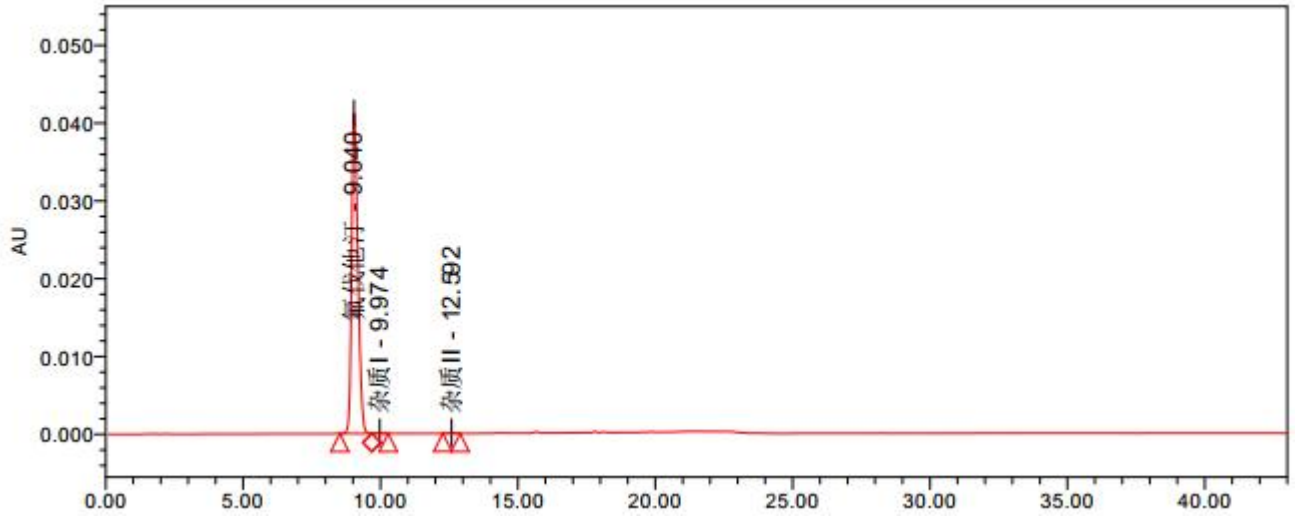
365nm

### 3. Conclusion (结论)

Using Welch Ultimate® Alk C18(4.6×150mm, 5μm), the detection requirements of the customer can be met under these chromatographic conditions.

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