Test report

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Sample Information 样品信息			
Sample Name 样品名称	Lepidii Semen 葶苈子	Sample properties 样品性状	Yellowish-brown powder 黄棕色粉末
Date of receipt 收样日期	2026/01/12	During testing 测试期间	2026/01/20~01/23
Test the ingredients and structural formula 测试成分及结构式			
Quercetin-3-O-β-D-glucose--7-O-β-D-gentian disaccharide glycosides 槲皮素-3-O-β-D-葡萄糖-7-O-β-D-龙胆双糖苷			
Experimental Requirements 实验要求			
The theoretical plate count calculated based on the peak of quercetin-3-O-β-D-glucose--7-O-β-D-gentian disaccharide glycosides should be no less than 5800. 理论板数按槲皮素-3-O-β-D-葡萄糖-7-O-β-D-龙胆双糖苷峰计算应不低于 5800。			
Reference Method 参考方法			
Pharmacopocia 药典			
Reagent information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Item Number 货号
Acetonitrile 乙腈	HPLC 色谱纯	Welch 月旭	00814-01010

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1. Test process:

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1.1. Chromatographic conditions:

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Chromatographic column: 色谱柱	Ultimate® ODS-3 (4.6×250mm, 5µm) 00275-31043
Mobile phase: 流动相	0.1% acetic acid / acetonitrile = 89/11 0.1%醋酸/乙腈=89/11
Flow rate: 流速	1.0mL/min
Injection volume: 进样量	25µL
Column temperature: 柱温	25°C
Detector: 检测器	UV
Detection wavelength: 检测波长:	254nm
Notes: 注意事项	<p>Under these chromatographic conditions, the sample contains relatively complex components, with strongly retained substances not fully eluted after approximately 100 minutes following the main peak. If conditions permit, it is recommended to add an additional gradient segment after the main peak to elute the strongly retained substances as much as possible, thereby minimizing their impact on subsequent injections</p> <p>该色谱条件下, 样品中成分较杂, 主峰后约 100 分钟还有强保留物质未被完全洗脱, 条件允许的情况下, 建议在主峰出峰后增设一段梯度, 将强保留物质尽可能洗脱下来, 减少对下一针的影响</p>

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制)

Take 1mL of acetic acid, add it to 1000mL of ultrapure water, mix well. Take 890mL of the above solution, add 110mL of chromatographic grade acetonitrile, shake well and degas by ultrasonic waves.

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1.2.2. Reference solution (对照品溶液)

Take an appropriate amount of reference substance, dissolve in 30% methanol and dilute to a solution of 20 μ g/mL.

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1.2.3. Test solution (供试品溶液):


Take 1.0g of the test sample powder, add 50mL of 70% methanol, heat and reflux for 1 hour, shake well and filter.

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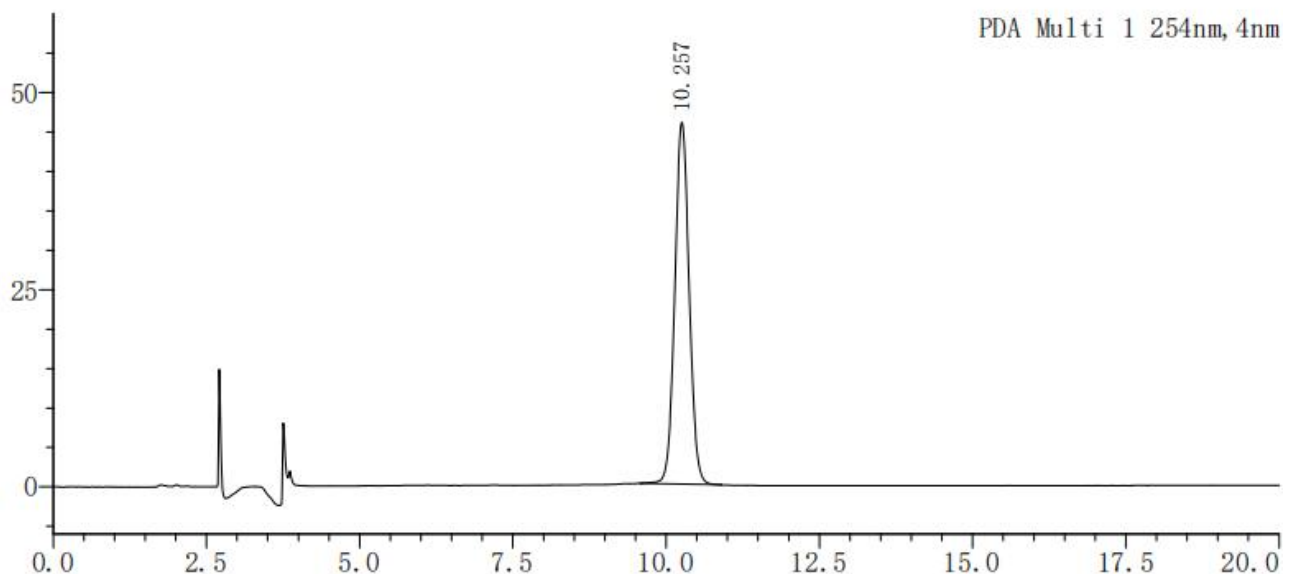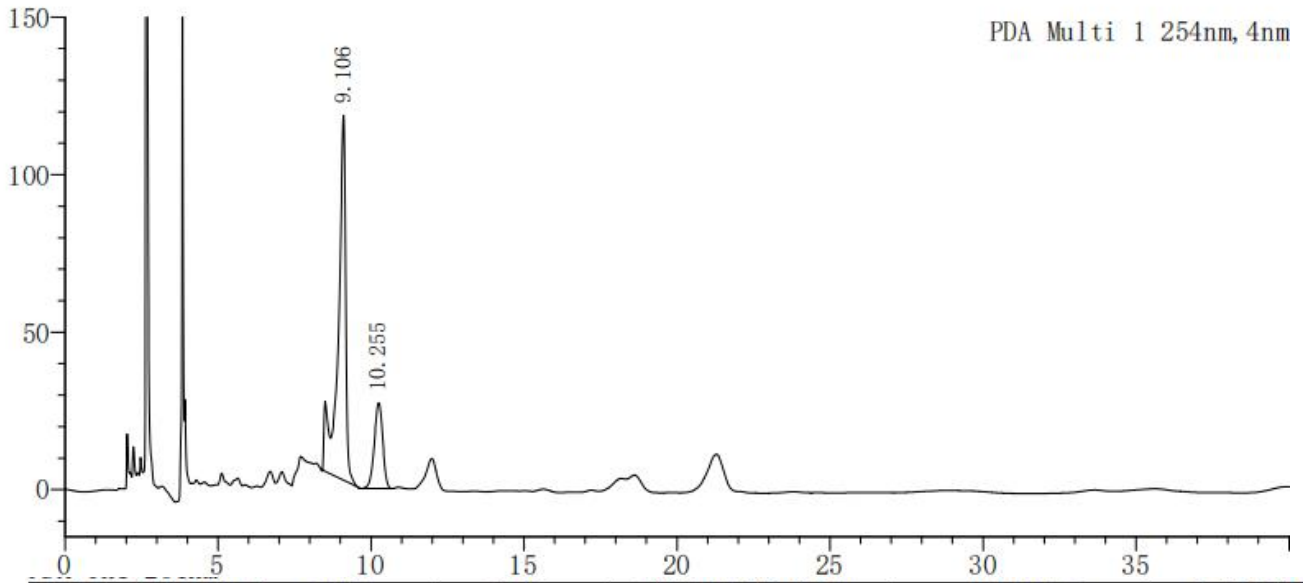
3. Conclusion (结论)

Using Welch Ultimate® ODS-3 (4.6×250mm, 5µm), the detection requirements of the customer can be met under these chromatographic conditions.

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