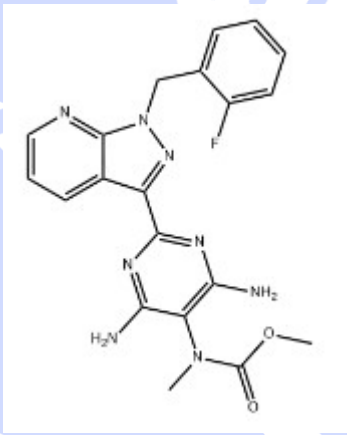
## Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Leo Siega 利奥西呱	Sample properties 样品性状	Solid 固体
Date of receipt 收样日期	2025/12/17	During testing 测试期间	2025/12/27 ~ 2025/12/29
Test the ingredients and structural formula			
测试成分及结构式			
 Leo Siega 利奥西呱			
Experimental Requirements			
实验要求			
The retention time of the main peak of the content reference solution should be stable, and the RSD of the peak area of the relevant reference solution after continuous injection of 6 needles should be less than 5.0%			
含量对照品溶液主峰保留时间需稳定, 有关对照品溶液连续进样 6 针峰面积 RSD 小于 5.0%			
Reference Method			
参考方法			
/			
Reagent information			

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## 1.2. Solution preparation (溶液配制)

### 1.2.1. Preparation of mobile phase (流动相配制):

Mobile Phase A: Weigh 2.3g of ammonium dihydrogen phosphate and dissolve it in 1000ml of water. Adjust the pH value to 3.5 with phosphoric acid. Measure 740ml, add 260ml of acetonitrile, mix well and filter to obtain the product.

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### 1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

Diluent: Phosphoric acid solution (pH1.6)- acetonitrile (80:20)

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### 1.2.3. Test solution (测试溶液):

1. Content reference substance solution: (Approximately 0.1mg of lioxigua per 1ml)

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2. Reference substance solution related to the substance: (Approximately 0.2µg of lioxigua per 1ml)

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3. System suitability solution: (Approximately 0.1mg of lioxigumab, 0.5µg/ml NK1 reference substance and 0.5µg/ml NK2 reference substance per 1ml)

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4. Sensitivity solution: (Approximately 0.1µg of lioxipa per 1ml)

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Web: www.welchmat.com

## 2. Chromatogram and data (谱图和数据)

Analysis spectrum of Test solution (测试溶液分析图谱):

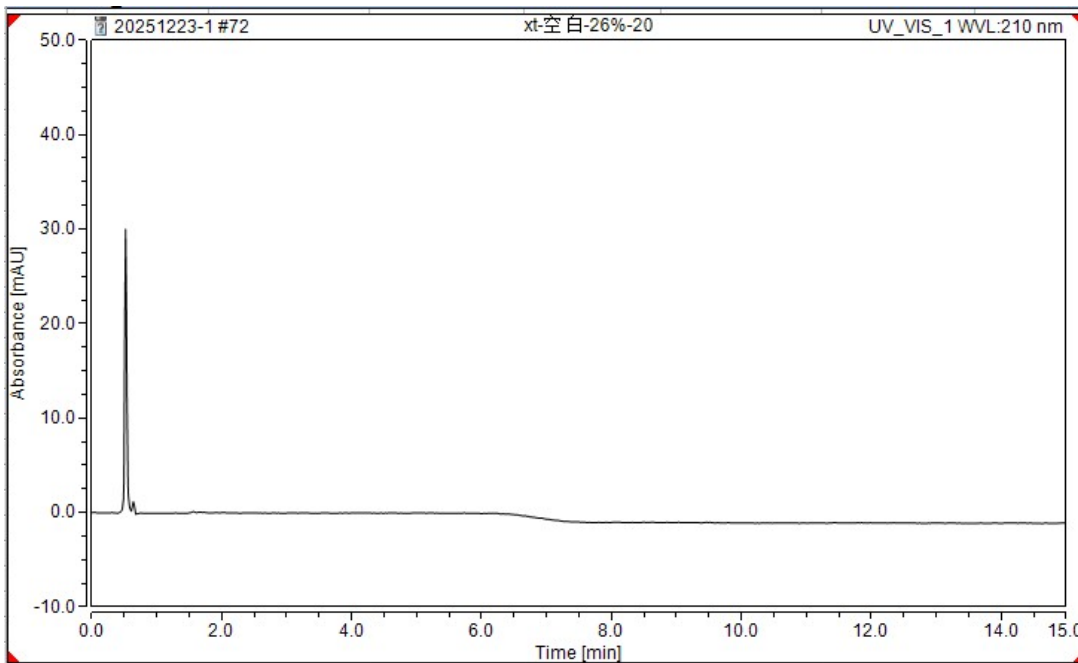
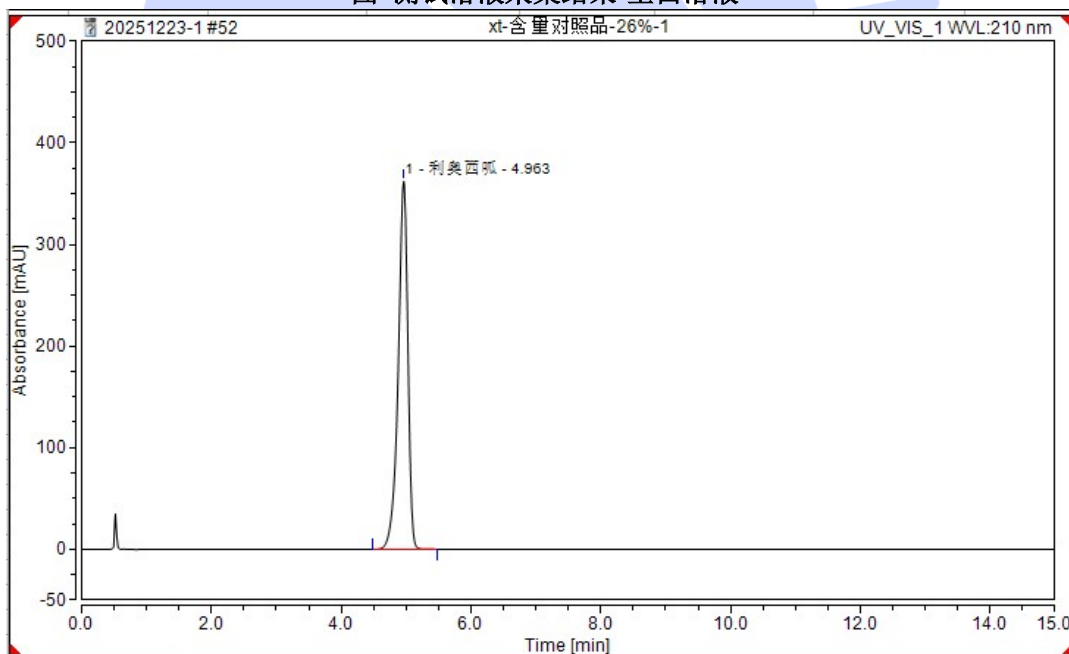
Figure : Acquired Results of the Test Solution-Blank solution  
图 测试溶液采集结果-空白溶液

Integration Results							
No.	Peak Name	Retention Time min	Area mAU*min	Height mAU	Resolution (EP)	Asymmetry (EP)	Plates (EP)
n.a.	NK1	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
1	利奥西呱	4.963	62.265	362.300	n.a.	0.82	5805
n.a.	NK2	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
<b>Total:</b>			<b>62.265</b>	<b>362.300</b>	<b>0.00</b>	<b>0.82</b>	

Figure : Acquired Results of the Test Solution-Content reference substance solution  
图 测试溶液采集结果-含量对照品溶液

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Web:www.welchmat.com

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Report No (报告编号): SHGZL-20251212-008

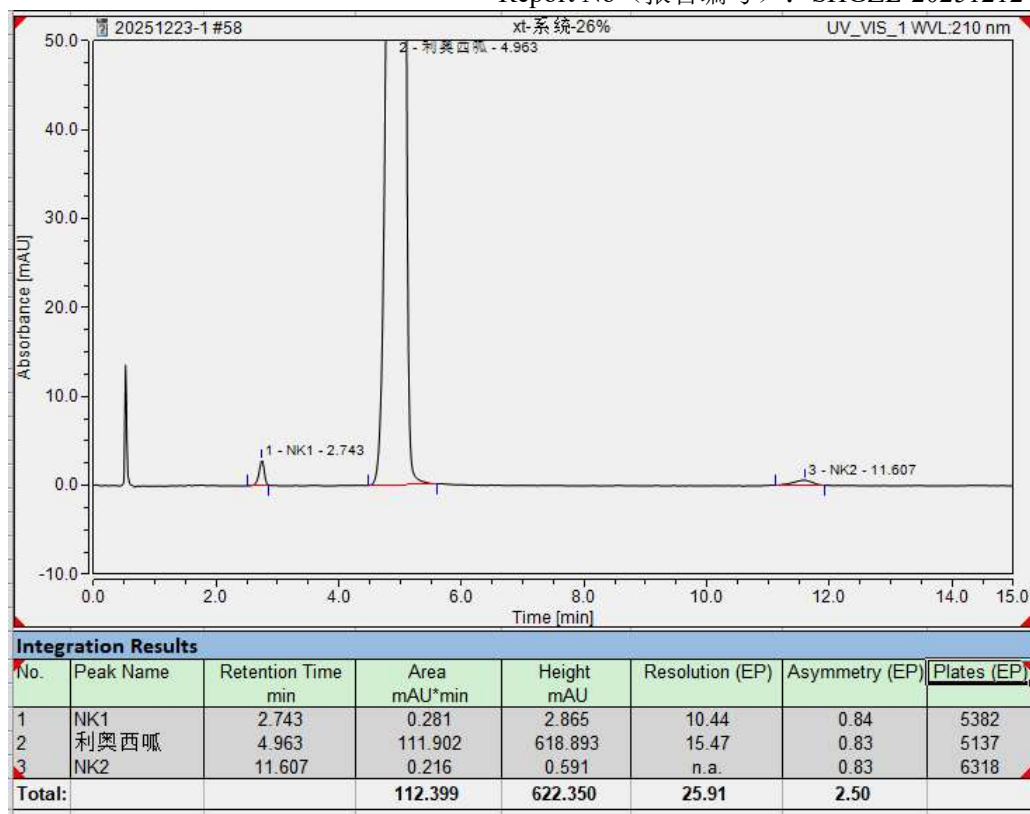
Figure : Acquired Results of the Test Solution-System suitability solution  
图 测试溶液采集结果-系统适用性溶液

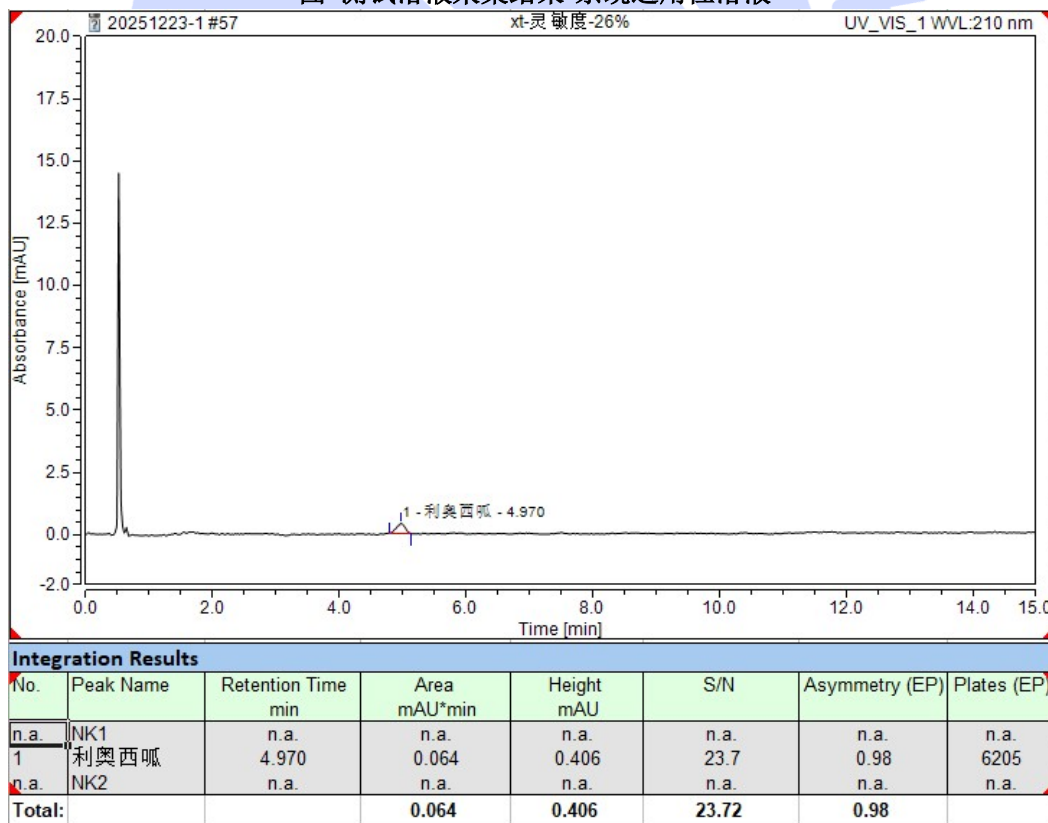
Figure : Acquired Results of the Test Solution-Sensitivity solution  
图 测试溶液采集结果-灵敏度溶液

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Tel:400-810-6969

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Web:www.welchmat.com

Test the analysis results of the solution (测试溶液分析结果)

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### 3. Conclusion (结论)

Using the Welch Materials Ultimate LP-C18 chromatographic column (4.6×250 mm, 5 μm) under these chromatographic conditions, The impurity separation degree before and after the main peak is both greater than 1.5, which can effectively separate the impurities before and after the main peak.

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