Test Report

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Sample Information 样品信息			
Sample Name 样品名称	Epimedii Folium 淫羊藿	Sample Trait 样品性状	brown solid powder 褐色固体粉末
Date of Receipt 收样日期	2026/01/05	Test Period 测试期间	2026/02/28~03/03
Target Analyte Structure 测试成分及结构式			
Icariin 淫羊藿苷			
Experimental Requirements 实验要求			
<p>Using icariin reference standard as the reference peak (S), the relative retention times of epimedin A, epimedin B, and epimedin C are approximately 0.73, 0.81, and 0.90, respectively, and should be within $\pm 5\%$ of the specified values.</p> <p>以淫羊藿苷对照品为参照 S 峰, 朝藿定 A、朝藿定 B、朝藿定 C 峰的相对保留时间分别为 0.73, 0.81, 0.90, 其相对保留时间应在规定值的$\pm 5\%$范围之内。</p>			
Reference Method 参考方法			
ChP 2025 Volume I 2025 版中国药典 一部			
Reagent Information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Item Number 货号

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1.2. Solution Preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制)

A: Take an appropriate amount of ultrapure water and degas by ultrasonication.

B: Take an appropriate amount of HPLC-grade acetonitrile and degas by ultrasonication.

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1.2.2. Standard solution (对照品溶液)

Take an appropriate amount of icariin reference standard, add methanol to prepare a solution containing 40 $\mu\text{g}/\text{mL}$.

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1.2.3. Test solution (供试品溶液)


Weigh about 0.2 g of the powder, place it in a stoppered conical flask, add 20 mL of dilute ethanol, sonicate for 1 h, mix well, filter, and collect the filtrate.

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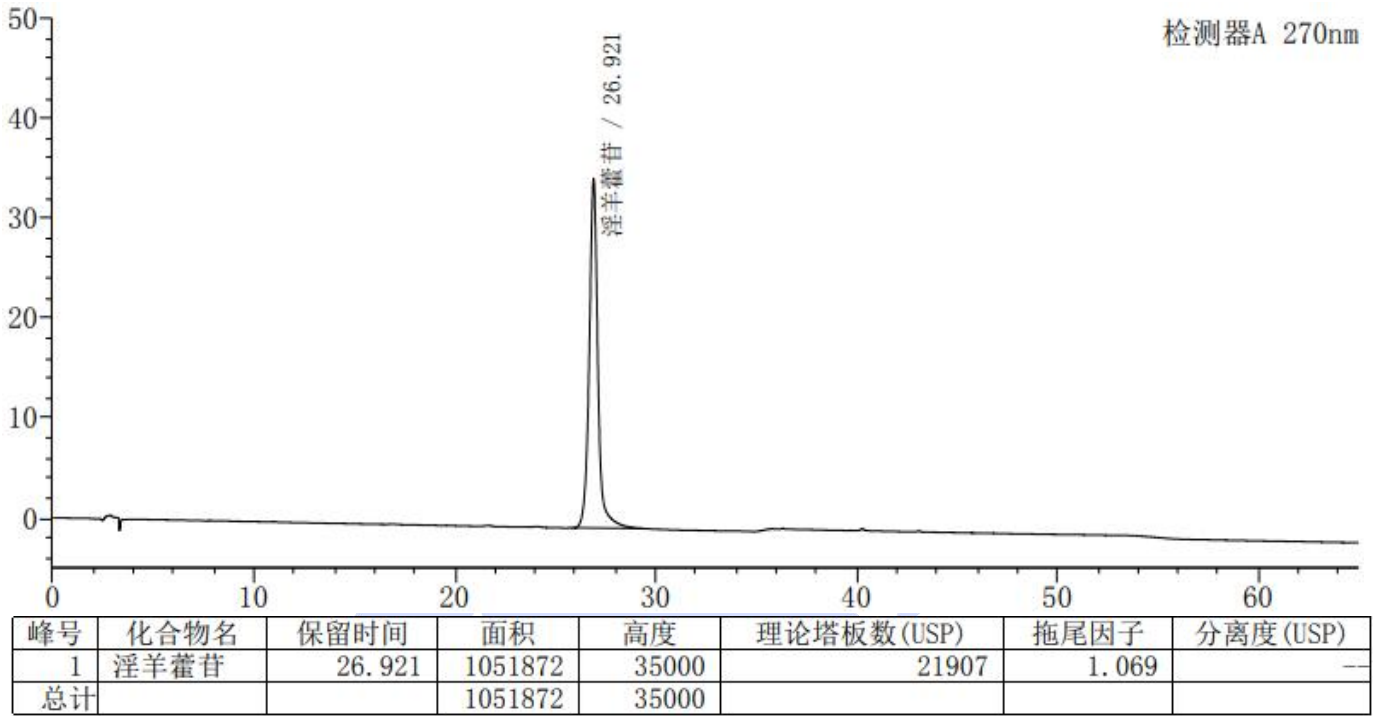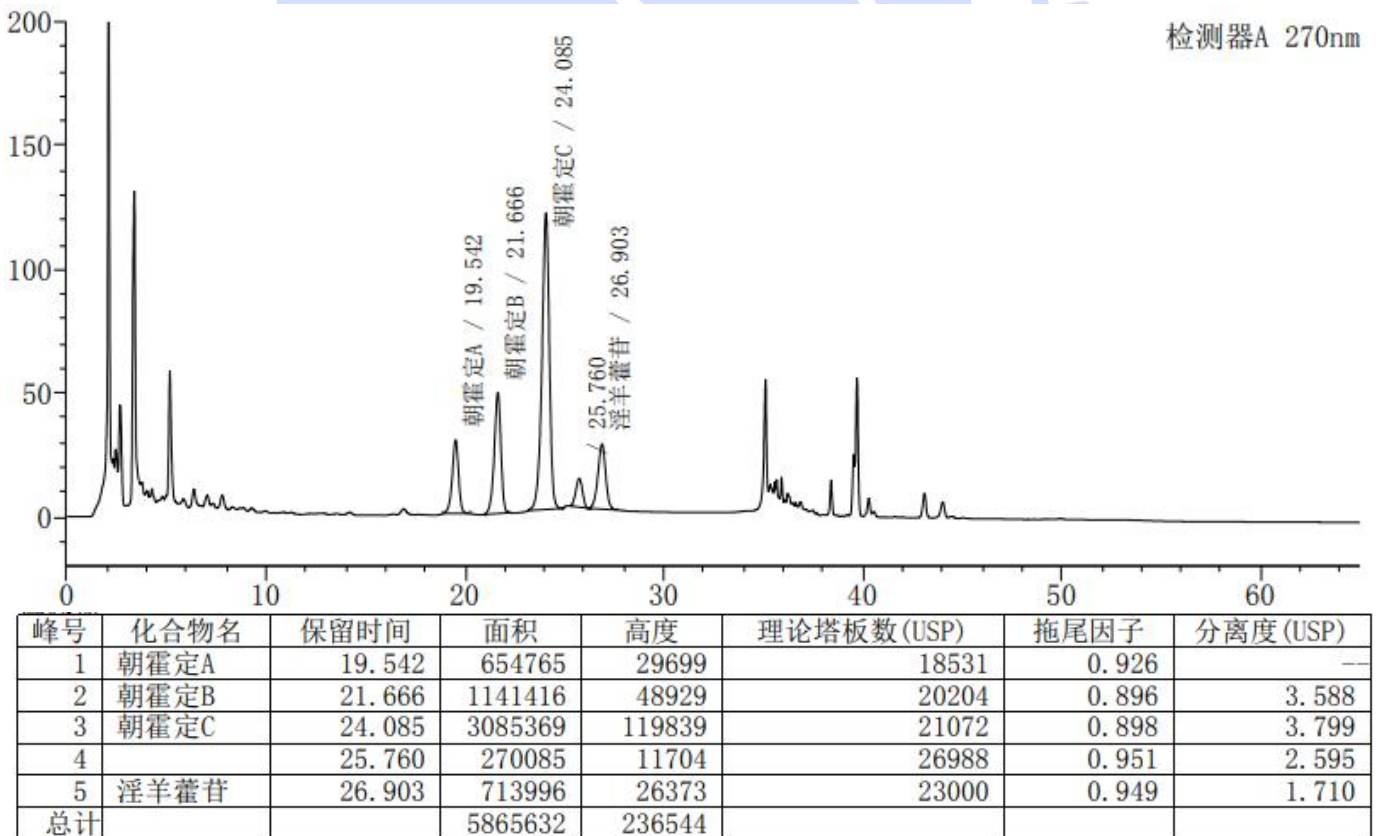
3. Conclusion (结论)

Using Welch Xtimate® C18 (4.6×250 mm, 5 μm), the detection requirements of the customer can be met under these chromatographic conditions.

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