Test report

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Sample Information 样品信息			
Sample Name 样品名称	Neostigmine Methylsulfate 甲硫酸新斯的明	Sample properties 样品性状	white solid powder 白色固体粉末
Date of receipt 收样日期	2025/12/19	During testing 测试期间	2025/12/24~01/05
Test the ingredients and structural formula 测试成分及结构式			
<p>氨基甲酸酯杂质/杂质 C[Ⓢ] 3-(二甲基)苯基二甲氨基甲酸酯 C₁₁H₁₆N₂O₂ 208.26[Ⓢ]</p> <p>羟基杂质/杂质 A[Ⓢ] 3-羟基-N,N,N',N'-三甲基苯胺甲硫酸盐 C₁₀H₁₁NO₃S 263.31[Ⓢ]</p> <p>杂质 D[Ⓢ] N,N'-二甲氨基甲酸3-(甲基氨基)苯酯 C₁₀H₁₄N₂O₂ 194.23[Ⓢ]</p> <p>苯酚杂质/杂质 B[Ⓢ] 3-(二甲胺基)苯酚 C₈H₁₁NO 137.18[Ⓢ]</p> <p>二甲氨基甲酸[Ⓢ]</p>			
Experimental Requirements 实验要求			
<p>Development method to ensure that the main components and impurities A, B, C, and D do not interfere with the peak of dimethylaminoformic acid, and the blank is interference free.</p> <p>开发方法, 使主成分以及杂质 A、B、C、D 不干扰二甲氨基甲酸出峰, 空白无干扰。</p>			
Reference Method 参考方法			
None 无			
Reagent information			

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制)

A: Take 1.36g of potassium dihydrogen phosphate, add it to 1000mL of ultrapure water, dissolve and filter, adjust the pH to 3.0 with phosphoric acid, and degas with ultrasound to obtain.

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B: Take an appropriate amount of chromatographically pure acetonitrile and sonicate to obtain it.

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60% acetonitrile water.

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1.2.3. Single standard solution (单标溶液)

Single standard solutions of main components and impurities A, B, C, and D: Take appropriate amounts of each single standard provided by the customer, and prepare solutions with impurity concentration of 1mg/mL and main component concentration of 3mg/mL using 60% acetonitrile water.

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Dimethyl aminoformic acid solution: Take an appropriate amount of the single standard provided by the customer and prepare a solution with a concentration of 5mg/mL using 60% acetonitrile water.

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1.2.4. Mixed standard solution (混标溶液)

Mixed standard solution 1: Take 0.5mL of the main component solution, 50uL of impurities A, B, C, and D each, and add 60% acetonitrile to 1mL to obtain.

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Mix standard 2 solution: Take 0.5mL of mixed standard 1 solution and 0.5mL of dimethyl aminoformic acid solution, mix well to obtain.

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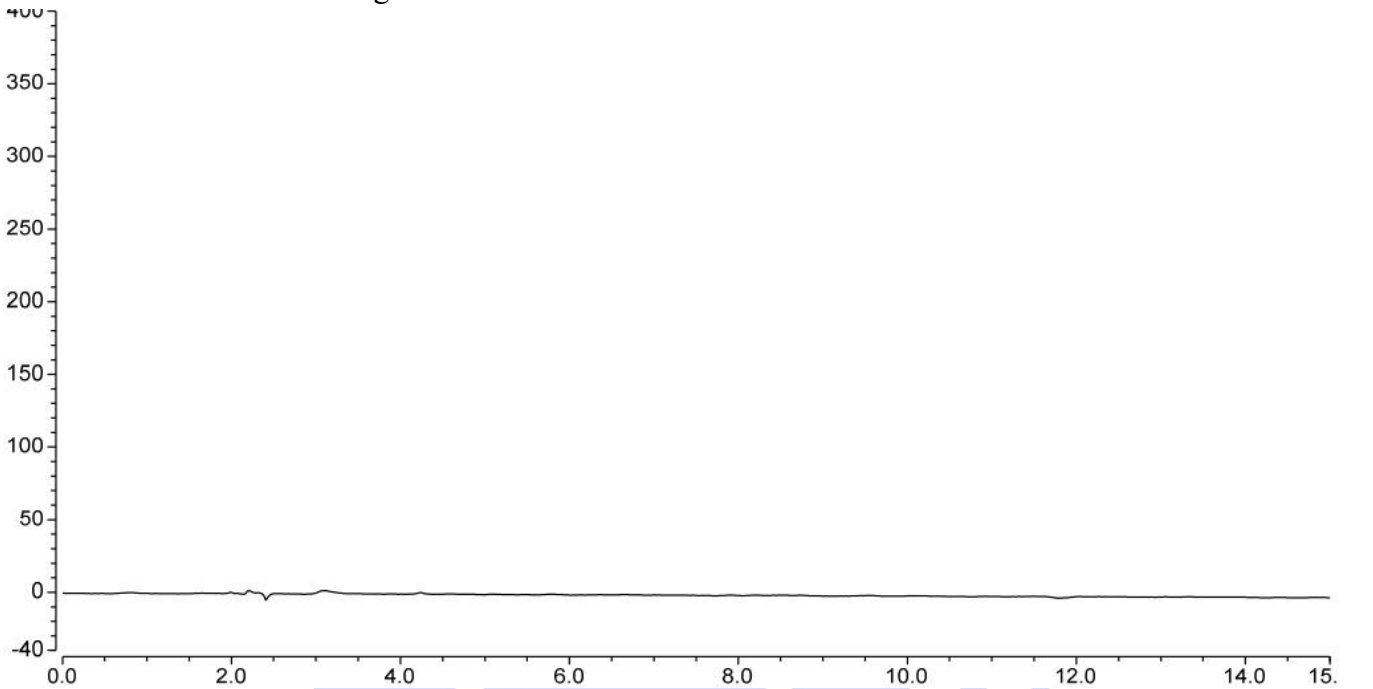
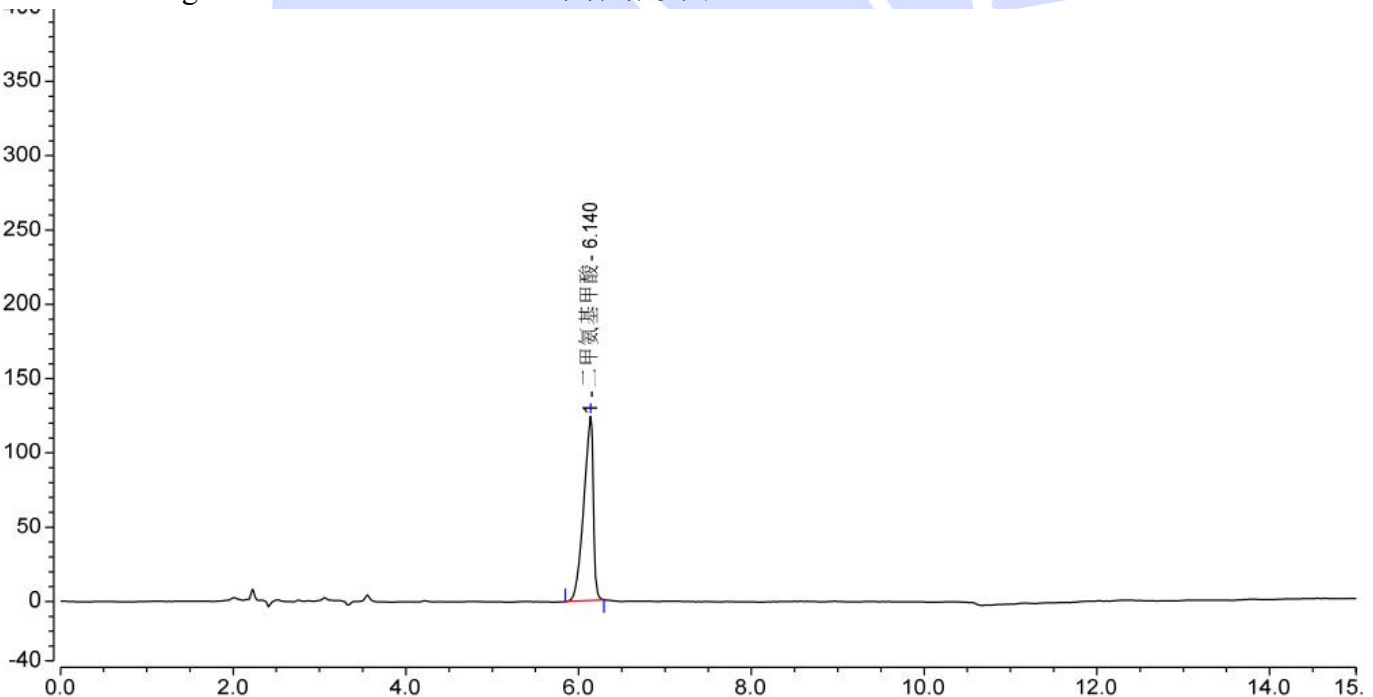
(2) Single standard solution view (单标溶液图)

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