Test report

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Sample Information			
样品信息			
Sample Name 样品名称	Pyrantel Pamoate 双羟萘酸噻嘧啶	Sample properties 样品性状	White solid powder 白色固体粉末
Date of receipt 收样日期	2026/01/14	During testing 测试期间	2026/01/19
Test the ingredients and structural formula			
测试成分及结构式			
Pyrantel Pamoate、Dihydroxynaphthalene acid 双羟萘酸噻嘧啶、双羟萘酸			
Experimental Requirements			
实验要求			
The method remains unchanged, and the retention time of dihydroxynaphthalene acid in the control sample is consistent with that of the salt forming dihydroxynaphthalene acid pyrimidine in the test sample. 方法不改动, 对照品中的双羟萘酸与供试品中成盐的双羟萘酸噻嘧啶的双羟萘酸保留时间一致			
Reference Method			
参考方法			
2025 Pharmacopoeia Volume 2 2025 版药典二部			
Reagent information			
试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	Item Number 货号
Acetonitrile	HPLC	Welch	00814-01010

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Web: www.welchmat.com

1.2.1. Preparation of mobile phase (流动相配制)

Take 30mL of glacial acetic acid and 12mL of diethylamine and add them to 928mL of chromatographically pure acetonitrile and 30mL of ultrapure water. Mix well and degas with ultrasound to obtain the product.

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1.2.2. Dihydroxynaphthalene acid solution (双羟萘酸溶液)

Take an appropriate amount of solid control provided by the customer and mix it with a flow mixture to form a solution of 0.08mg/mL.

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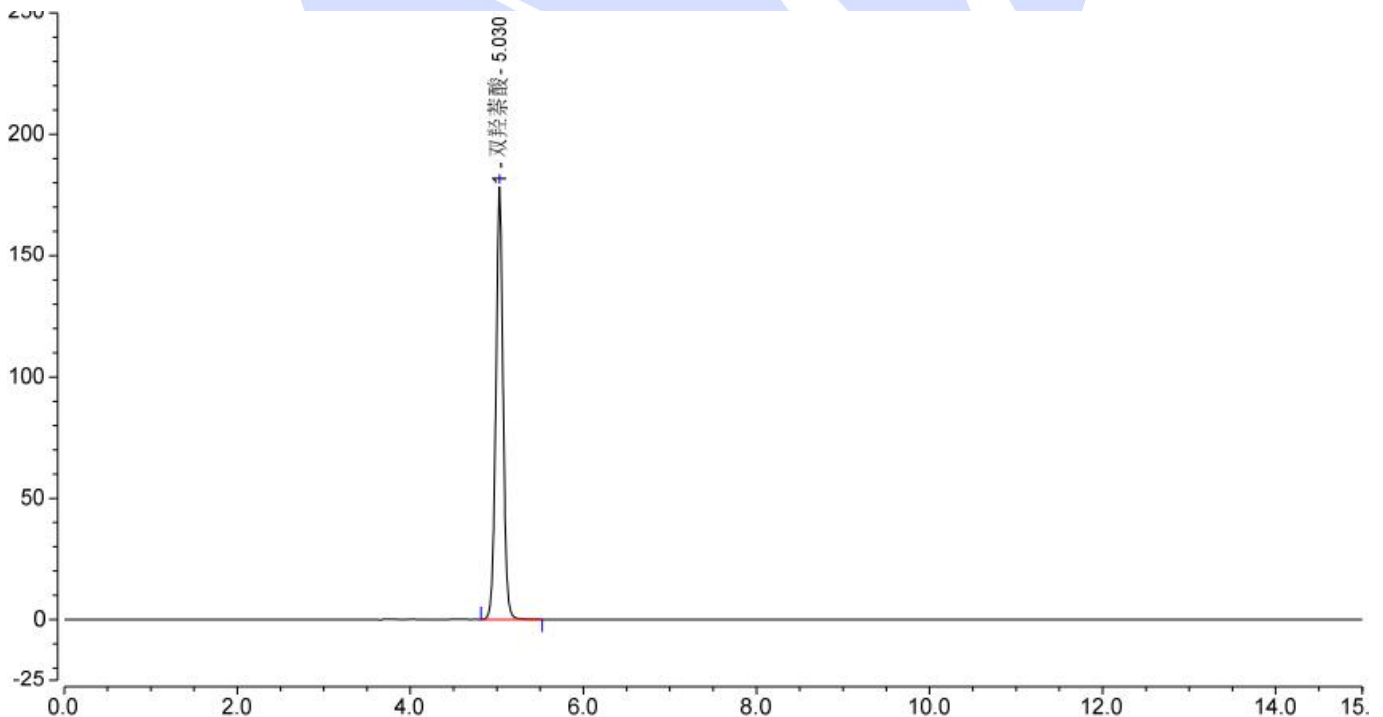
1.2.3. Thiopyrimidine dihydroxynaphthalene solution (双羟萘酸噻嘧啶溶液)

Take an appropriate amount of solid control provided by the customer and mix it with a flow mixture to form a solution of 0.08mg/mL.

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2. Chromatography and data (谱图和数据)

(1) Dihydroxynaphthalene acid solution view (双羟萘酸溶液图)

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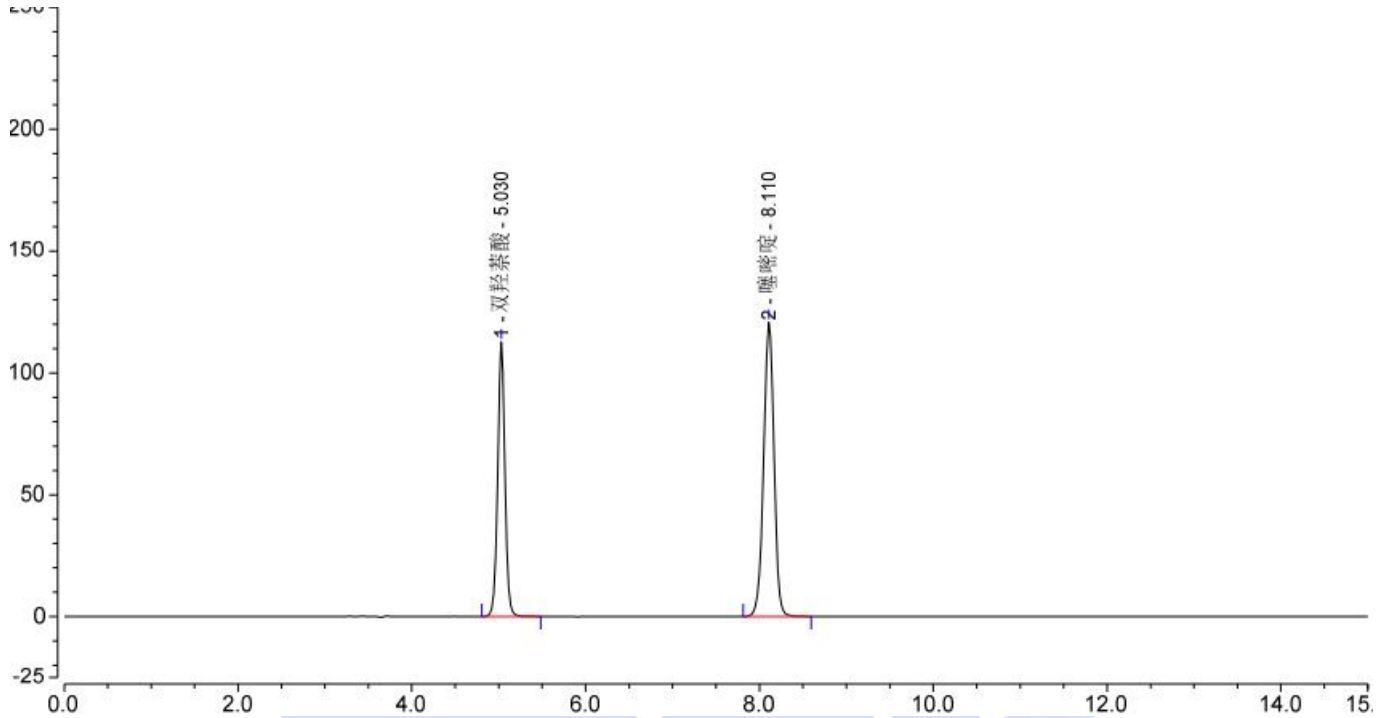
(2) Thiopyrimidine dihydroxynaphthalene solution view (双羟萘酸噻嘧啶溶液图)

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3. Conclusion (结论)

Using Welch Ultimate® HILIC-Silica (4.6×250mm,5µm), the detection requirements can be met under these chromatographic conditions.

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